

ChemiDoc and ChemiDoc MP Imaging Systems with Image Lab Touch Software

User Guide

Version 2.4



BIO-RAD

ChemiDoc and ChemiDoc MP Imaging Systems with Image Lab Touch Software

User Guide

Version 2.4

The Bio-Rad logo is located at the bottom center of the page. It consists of the words "BIO-RAD" in a bold, white, sans-serif font, set against a dark, rounded rectangular background.

BIO-RAD

Technical Support Department

The Technical Support department in the U.S. is open Monday through Friday, 5:00 AM to 5:00 PM, Pacific time.

Phone: 1-800-424-6723, option 2

Email: Support@bio-rad.com (U.S./Canada Only)

For technical assistance outside the U.S. and Canada, contact your local technical support office or click the Contact us link at www.bio-rad.com.

Notice

No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopy, recording, or any information storage or retrieval system, without permission in writing from Bio-Rad.

Bio-Rad reserves the right to modify its products and services at any time. This guide is subject to change without notice. Although prepared to ensure accuracy, Bio-Rad assumes no liability for errors or omissions, or for any damage resulting from the application or use of this information.

Image Lab Touch software is based in part on the work of the CImg project (<http://cimg.eu/>). See license for details at www.cecill.info/licences/Licence_CeCILL-C_V1-en.html.

Image Lab Touch is based in part on libraries from GCC runtime and the Gnu C library.

ChemiDoc and ChemiDoc MP Imaging Systems use the Linux operating system, which contains software licensed under the following licenses as well as others:

- Gnu Public License v 1.0, 2.0, 3.0
- Lesser Gnu Public License v 2.0, 2.1, 3.0
- Apache License 2.0
- BSD License
- Artistic

Clarity Max Western ECL Substrate is manufactured by Cyanagen Srl and is the subject of patent application numbers US7855287, EP1950207, US9040252, AU2011202658, CA2742025, US8129136, and EP1962095, together with other equivalent granted patents and patent applications in other countries like CN102313732.

SYBR is a trademark of Thermo Fisher Scientific Inc. Bio-Rad Laboratories, Inc. is licensed by Thermo Fisher Scientific Inc to sell reagents containing SYBR Green I for use in real-time PCR, for research purposes only.

Bio-Rad, Mini-PROTEAN, and Trans-Blot are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions.

All trademarks used herein are the property of their respective owner.

Copyright © 2019 by Bio-Rad Laboratories, Inc. All rights reserved.

Table of Contents

Safety and Regulatory Compliance	11
Important Safety Information	11
Warranty	12
General Precautions	13
Regulatory Notices	13
Safety Alerts	14
Types of Safety Hazards	14
Levels of Potential Risk	15
Instrument Safety Warnings	16
Notice	17
Power Safety	17
Fuses	17
Chapter 1 Introduction	19
Product Features	20
Front Panel Components	21
Rear Panel Components	22
CCD Camera	23
Image Lab Touch Software	23
Emission Filters	23
Optional Accessories	23
Printer	23

Sample Trays	23
Technical Specifications	25
Environmental Requirements	28
Chapter 2 Starting Image Lab Touch	29
About the Welcome Screen	30
Interacting with the Touch Screen	31
Entering Text on the Touch Screen	32
Choosing Imaging Settings	34
Locking the System	36
Unlocking the System	37
Getting Help	38
Chapter 3 Logging In to Image Lab Touch	39
Logging in as a User	39
Creating a Password	40
Removing a Password	43
Logging in as an Administrator	43
Setting Administrator Options	46
Chapter 4 Understanding the Views	49
Main Toolbar Settings	49
Camera View	51
Preview	53
Image View	54
The Image View Toolbar	55
The Gallery	56
Color Cues	58

Chapter 5 Choosing a Sample Tray	63
Applications Listed by Tray Type	65
Blot/UV/Stain-Free Tray	65
Blue Tray	68
White Tray	68
Chapter 6 Acquiring an Image	73
Placing the Sample in the Imager	74
Acquiring an Image — General Steps	74
Setting Image Type and Size	76
Choosing Image Type	76
Choosing Image Size	76
Choosing the Application	78
(Multichannel) Selecting a Channel Color	80
Previewing an Image	81
Exposure Settings	83
Exposure Settings for Most Gels and Blots	83
Exposure Settings for Stain-Free Gels	84
Setting Exposure Time for a Chemiluminescent or Low-Light Blot	85
Automatic Exposure Settings	86
Setting Exposure Automatically for Chemiluminescent or Low-Light Applications	86
Setting the Exposure Manually	90
(Single Channel) Acquiring Signal Accumulation Mode (SAM) Images	92
Adjusting SAM Acquisition Settings	95
Acquiring the Image	96
(Stain-Free Gel) Skipping Gel Activation	97
(Stain-Free Gel) Acquiring an Image after Canceling Acquisition	98
(Multichannel) Retaking an Image	98

Merging Images	101
Naming Merged Images	103
Chapter 7 Excising Bands from Samples	105
Excising Bands on a White or Blue Tray	105
Excising Bands on a Blot/UV/Stain-Free Tray	106
Chapter 8 Viewing Images	111
Accessing the Gallery	111
Selecting and Opening Images in the Gallery	113
Filtering Gallery Images by Date	115
Checking Intensity Values	117
Viewing Multichannel Images in Image View	120
Viewing a Composite or Merged Image in Gray Scale	122
Comparing Images	123
Copying Image Settings	125
Viewing and Editing Image Information	126
Viewing SAM Images	128
Chapter 9 Managing Images	131
Adding Notes to an Image File	131
Changing the File Naming Format	132
Renaming Images	134
Renaming SAM Images	135
Printing an Image	136
Searching for Images in the Gallery	137
Deleting Images	137
Deleting SAM Images	138

Chapter 10 Exporting Images	141
Preparing to Export Images	142
Creating a Shared Folder on a Windows Computer	142
Creating a Shared Folder on a Mac Computer	145
Determining the UNC of a Previously Shared Folder	149
Export Formats	152
Exporting Images	155
Exporting to a USB Flash Drive or External Hard Drive	158
Exporting to a Shared Folder for the First Time	159
Exporting to a Shared Folder on a Network or Computer	161
Entering Log-in Credentials	161
Exporting to a Recent Location	163
Removing Network Locations	164
Disconnecting from a Shared Folder	164
Chapter 11 System Settings	165
Setting the System Date and Time	165
Setting the Sound Volume	167
Appendix A Maintaining the Imaging System	169
Reinstalling Image Lab Touch Software	169
Maintaining Transilluminator Bulbs	170
Replacing UV Transilluminator Bulbs	170
Replacing the Fuses	173
Appendix B Important Imaging Concepts	177
Exposure Settings for Chemiluminescent and Low-Light Applications	177
Adjusting How Images Are Displayed	178
Adjusting Image Brightness and Contrast	178

Table of Contents

Restoring Default Settings	181
The Image Resolution/Time Scale	182
How Binning Settings Affect Image Quality	183
Appendix C Ordering Information	187

Safety and Regulatory Compliance

Note: This equipment has been tested and found to comply with the limits for a Class A digital device pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radiofrequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment is intended for laboratory use only and if used in a residential area is likely to cause harmful interference in which case users will be required to correct the interference at their own expense.

Important Safety Information

Please read these instructions before operating the ChemiDoc MP or ChemiDoc imaging system.

These instruments are suitable for research use only. Therefore, they must be used only by specialized personnel who know the health risks associated with the reagents that are normally used with these instruments.



WARNING: The imaging of some applications involves UV illumination.

These instruments should be used only by trained personnel who know the health risks associated with the UV radiation normally associated with these instruments. Users should be trained on the appropriate personal protective equipment for working with UV light to minimize UV exposure.

To perform band excision using the blot/UV/stain-free tray, the transilluminator drawer is pulled out with the UV source enabled. This exposes the user to UV radiation, which can cause permanent damage to the eyes and skin. In its lowered position, the instrument's optional acrylic shield provides some UV protection. However, in its raised position, it does not provide complete protection to the user, and it does not protect others who are standing in the area around the imager.

To protect users and bystanders, these procedures must be followed:

- Protect all skin surfaces (including the neck, ears, and hands). Before performing band excision, the user and anyone near the imager must put on personal protective equipment including UV protective safety glasses, a face shield, lab coat, and gloves. A typical and reasonable expectation of use is three operations per user a day for three minutes each.
- Bystanders without protective gear must stand at least 1.5 meters (five feet) away from the imager and limit their exposure to no more than one hour per day.

Note: There is no exposure to UV radiation with the blue or white trays. No protective gear is necessary when excising bands with these trays.

Warranty

The ChemiDoc MP and ChemiDoc imaging systems are warranted against defects in materials and workmanship for one year. If any defect occurs in the instrument during this warranty period, Bio-Rad Laboratories, Inc. will repair or replace the defective parts at its discretion without charge. The following defects, however, are specifically excluded:

- Defects caused by improper operation
- Repair or modification done by anyone other than Bio-Rad Laboratories, Inc. or the company's authorized agent
- Use of spare parts supplied by anyone other than Bio-Rad Laboratories, Inc.
- Damage caused by accident or misuse

- Damage caused by disaster
- Corrosion caused by improper solvents or samples

General Precautions

- Read the user guide carefully.
- Use the instrument only for the intended purpose of gel and blot image acquisition in research laboratories.
- Connect the instrument to a grounded power source and to a circuit breaker.
- Do not pour liquids on or inside the instrument.
- Clean the sample tray after use.

Regulatory Notices

The ChemiDoc and ChemiDoc MP imaging systems are designed and certified to meet EN 61010, the internationally accepted electrical safety standard, EMC regulations, and TUV requirements. Certified products are safe to use when operated in accordance with this user guide. Do not modify or alter these instruments in any way. Modification or alteration of these instruments will

- Void the manufacturer's warranty
- Void the regulatory certifications
- Create a potential safety hazard







Bio-Rad Laboratories, Inc. is not responsible for any injury or damage caused by use of these instruments for purposes other than those for which they are intended or by modifications of the instruments not performed by Bio-Rad Laboratories, Inc., or an authorized agent.

Safety Alerts

Alert icons appear in cautionary and warning paragraphs in this guide to call attention to safety hazards.

Types of Safety Hazards

Most alert icons depict the relevant type of safety hazard.

Alert Icon	Explanation
	<p>General</p> <p>Indicates a potential hazard requiring special attention. This icon appears when the hazard or condition is of a general nature.</p>
	<p>Electrical Hazard</p> <p>Indicates a potential hazard requiring special attention when you are working with electricity or electrical equipment.</p>
	<p>Extreme heat and flammable materials</p> <p>Indicates a potential hazard requiring special attention when you are working with extreme heat and flammable materials.</p>
	<p>Radiation hazard</p> <p>Indicates a potential hazard requiring special attention when you are working with UV radiation.</p>

Levels of Potential Risk

Each alert icon appears in a paragraph type in this guide that indicates the level of potential risk posed by the hazard or action described.

Cautions

A caution alerts you to take or avoid a specific action that could result in loss of data or damage to the instrument. A caution can also indicate that minor or moderate injury might occur if the precaution against a potential hazard is not taken.

Example



Caution: With the exception of cleaning or replacing light bulbs, refer all servicing to qualified Bio-Rad personnel or their agents.

Warnings

A warning precedes an action that, if not followed correctly, could cause serious injury or death to the operator, serious or total loss of data, or serious damage to the instrument.

Example






WARNING! Radiation Hazard

Keep the UV shield open for as little time as possible.

Instrument Safety Warnings

Before you operate the instrument, carefully read the explanation for each safety icon.

Safety Icon	Explanation
	Caution: With the exception of cleaning or replacing light bulbs, refer all servicing to qualified Bio-Rad personnel or their agents. If you experience technical difficulties with the instrument, contact Bio-Rad to schedule service. The instrument should not be modified or altered in any way. Alteration voids the manufacturer's warranty and might create a potential safety hazard for the user.
	WARNING! If any interlock is defeated, there is a possibility of UV-B radiation hazard due to UV-B light exposure. Exercise caution when servicing the instrument.
	WARNING! This instrument must be connected to an appropriate AC voltage outlet that is properly grounded.

Notice

Bio-Rad's imaging systems are intended for laboratory use only. These devices are meant for use by specialized personnel who know the health risks associated with reagents used in electrophoresis. The UV light source is computer controlled, and proper interlocks are implemented to avoid users' accidental exposure to UV radiation. Bio-Rad Laboratories, Inc. is not responsible for any injury or damage caused by use of these instruments for purposes other than those for which they are intended, or for instrument modifications not performed by Bio-Rad Laboratories, Inc. or an authorized agent.

Power Safety

The imaging systems use a universal power supply that automatically chooses the correct voltage for your country or region.

The operating voltage requirement for the ChemiDoc and ChemiDoc MP imaging systems is 100–240 VAC; 50–60 Hz.

Fuses

The imaging systems have two user-serviceable fuses, F1 and F2, which are located on the rear panel and are a part of the power entry module. See [Replacing the Fuses on page 173](#) for more information.

Chapter 1 Introduction

ChemiDoc and ChemiDoc MP imaging systems facilitate high sensitivity and high quality image acquisition for selected applications (gels and blots). Both systems use Bio-Rad's Image Lab Touch software to control image capture and optimization. You interact with the instrument via an integrated touch screen and a simple user interface. Using Image Lab Touch, you can acquire and view images, fine tune their appearance, and print them. You can also export images to a computer running Image Lab desktop software, in which you can analyze the images in detail.

Note: Image Lab Touch software does not support image analysis. Use Image Lab desktop software version 5.2 or greater on a separate computer to analyze images acquired with the imaging systems.

The Bio-Rad imaging systems are compact gel/blot imaging instruments. These instruments automate the process of selecting blot detection parameters and acquire high-quality and high-sensitivity gel and multichannel western blot images with the tap of an on-screen button.

- Using the ChemiDoc MP imaging system, you can acquire data from 1–3 channels in a multichannel image for applications that use blots, including fluorescence, chemiluminescence, and colorimetric applications.
- Using the ChemiDoc imaging system, you can acquire a two-channel image consisting of data from a chemiluminescent channel and a colorimetric channel, in addition to acquiring single-channel images.

Note: This user guide includes examples of images acquired with applications available only on ChemiDoc MP imaging systems. The accompanying instructions are accurate for both systems.


Product Features

The Bio-Rad imaging systems support the acquisition of images from a wide range of laboratory gels and blots. The instruments use a supersensitive camera with a charge-coupled device (CCD) and a large maximum aperture lens that provides highest chemiluminescent sensitivity. The imagers have a built-in UV transilluminator and white LEDs for epi (reflective) illumination. The ChemiDoc MP includes five additional LEDs for detection of a larger range of fluorophores and dyes. The imagers work with gels and blots stained with a wide range of dyes and fluorophores.

Additional features include


- Support for colorimetric western blotting applications, chemiluminescent imaging, and DNA/protein gel visualization
- Multiplex fluorescent western blotting applications (MP only)
- Blue, green, red, and infra red excitation sources for multiplex western blotting (MP only)
- Dynamic flat fielding specific to each application
- Smart, tray-based imaging that identifies the correct applications and presents appropriate filter and illumination sources for each
- Dynamic range of >4 orders of magnitude

Front Panel Components

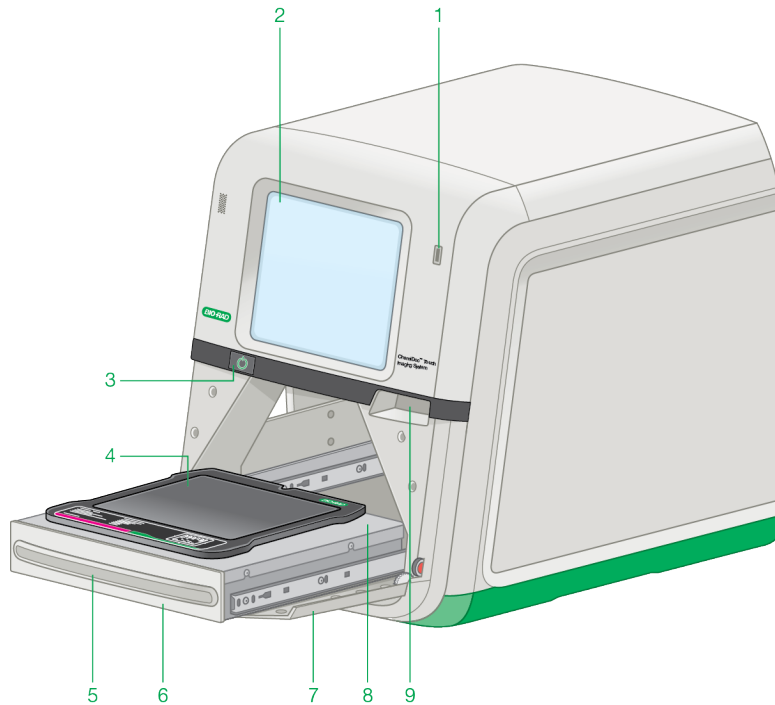

1	USB port	6	Transilluminator drawer
2	Touch screen	7	Front door (open position)
3	System on/off button	8	Imaging stage
4	Sample tray	9	Front door open handle
5	Transilluminator drawer handle		

Rear Panel Components

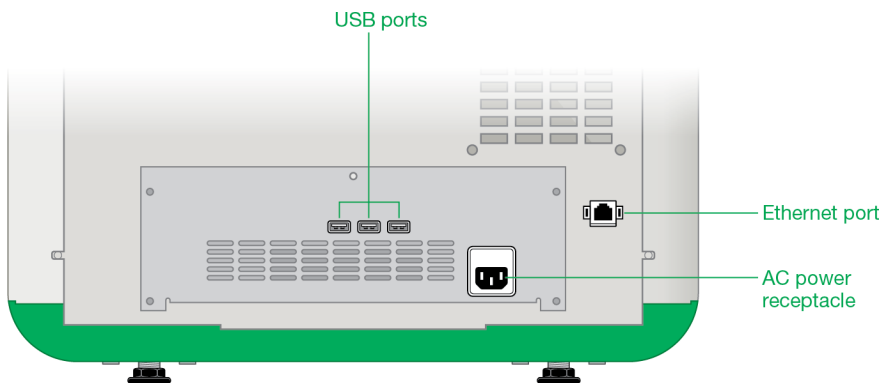

USB Ports

Use these ports to connect any of the following accessories to the imaging system.

- Printer
- Mouse
- Keyboard
- USB drive

Ethernet Port

Use this port to transfer files to a network drive.

Note: The ChemiDoc imaging system cannot be controlled by an external computer.

CCD Camera

Based on the application and imaging area you select, Image Lab Touch uses a patented algorithm to adjust the system camera/lens settings (iris, focus) automatically. See [Technical Specifications on page 25](#) for more information.

Image Lab Touch Software

The imager ships with Image Lab Touch software installed.

Emission Filters

The imagers have a motorized emission filter wheel that automatically moves to the correct position for the chosen application.

Optional Accessories

See [Appendix C, Ordering Information](#), for a list of optional accessories and replacement parts.

Printer

An optional USB printer, the Mitsubishi P95 thermal printer (cat. #1708089), is available from Bio-Rad for use with the imaging system.

Sample Trays

Blot/UV/Stain-Free Sample Tray

The blot/UV/stain-free tray (cat. #12003028) is shipped with the imaging systems. The UV tray is made of UV transmissive glass that blocks other wavelengths of light from the

transilluminator fluorescent bulbs, reducing the background light in the emission spectrum of samples imaged using trans-UV excitation. The blot/UV/stain-free tray can be used with all blotting applications.

White Sample Tray

The optional white tray (cat. #12003026) is for use with transillumination of colorimetric gels, such as coomassie or silver stained samples.

Blue SampleTray

The optional blue tray (cat. #12003027) is a UV-to-blue-light conversion screen that makes appropriately stained DNA samples visible while protecting them from UV damage.

See [Chapter 5, Choosing a Sample Tray](#), for more information.

Technical Specifications

Supported Applications	
Fluorescence	ChemiDoc MP only
Chemiluminescence	Yes
Colorimetry	Yes
Gel documentation	Yes
Hardware Specifications	
Touch screen functionality	<ul style="list-style-type: none">■ Multitouch capable■ Display resolution 1024 x 768 pixels■ 12.1" (30.73 cm) display
Onboard computer system	<ul style="list-style-type: none">■ 4 GB RAM■ 120 GB disk space■ 4 USB ports
Sample thickness	Maximum supported thickness: 5 mm
Maximum image area	<ul style="list-style-type: none">■ Length: 16.8 cm■ Width: 21 cm

Excitation source	<ul style="list-style-type: none">■ Trans-UV 302 nm (standard)■ Epi-white (standard)■ Trans-white (optional)■ Trans-blue (optional) LEDs for multiplex fluorescence detection (optional, MP only) <ul style="list-style-type: none">■ Blue■ Green■ Red■ Far Red■ Infra Red
Detector	Deeply cooled CCD
Pixel size	4.54 x 4.54 μm
Cooling system	Thermoelectric
Filter wheel	Automated
Emission filters	535–645 nm (standard) Additional emission filters for MP only <ul style="list-style-type: none">■ Chemi■ 532/28■ 595/36■ 715/30■ 835/50
Dynamic range	>4.0 orders of magnitude
Pixel density (gray levels)	65,535

Instrument size	<div><div></div>Depth: 62.6 cm (25")<div></div>Width: 50 cm (20")<div></div>Height: 54.3 cm (21")</div>
Instrument weight	32 kg (78 lb)
Operating Ranges	
Operating voltage	100–240 VAC, 50–60 Hz
Operating temperature	10–28°C
Operating humidity	10–80% relative humidity (noncondensing)
Automation Capabilities	
Workflow automated selection	Application driven, tray-based imaging Application-driven autoselection of emission filter and excitation source
Autofocus	Precalibrated focus for any zoom setting
Image flat fielding	Dynamic; precalibrated and optimized per application
Autoexposure	2 user-defined modes (rapid or optimal)

Environmental Requirements

The imaging systems require a space 52 x 56 x 62 cm (W x H x D) and a clearance of at least 8 cm from the back for instrument ventilation and for connecting or disconnecting the AC power cord. Place the imager on a sturdy and level laboratory bench or table away from excessive heat and moisture. The imager's operating temperature range is 10–28°C. The imager contains a universal power supply that supports a voltage range of 100–240 VAC.




WARNING! Transilluminators are powerful sources of UV radiation, which can cause serious damage to unprotected eyes and skin. The accessory UV shield (cat. #1708375) provides UV protection. However, this shield does not guarantee complete protection nor does it provide protection to others in the area around the imager. Before performing band excision, the user and anyone near the imager must put on protective gear including eyeglasses (laboratory glasses provide adequate protection), a face shield, lab coat, and gloves.

Chapter 2 Starting Image Lab Touch

Press the On button on the imaging system. The imaging system turns on and Image Lab Touch software starts automatically.

The Welcome screen appears.

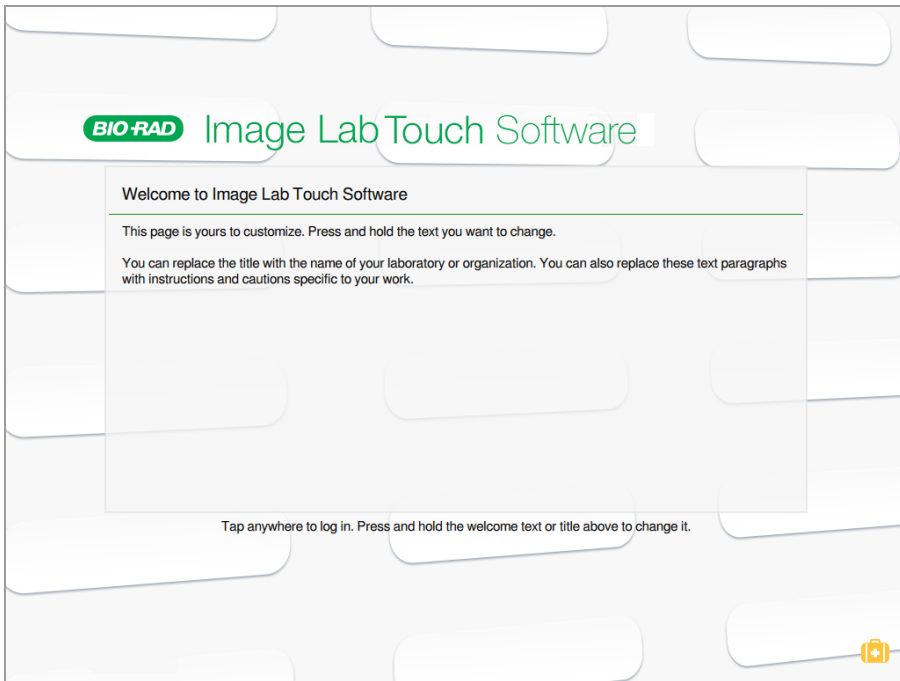

About the Welcome Screen

On the Image Lab Touch Software Welcome screen, you can customize the text for your organization. For example, you can enter the name of your business or laboratory in the upper box and add a warning message or other helpful information in the second box.

You can enter one line of text in the upper box. You can fill the lower box with text. The box does not scroll.

To edit Welcome screen text


1. Press and hold on the text you want to change.

The on-screen keyboard appears.

2. Enter your text.
3. Tap OK to save your changes.

What's in the Suitcase?

Tapping the Suitcase in the lower right corner of the Welcome screen displays two commands:

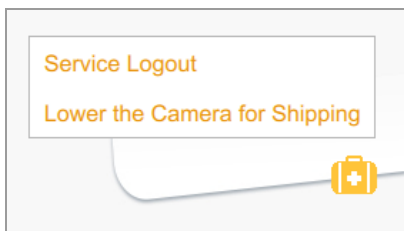

Service Logout — enables service personnel to log out of Image Lab Touch and access service tools.

Lower the Camera for Shipping — accesses detailed instructions for preparing the camera for shipment. For more information, see Repacking the ChemiDoc and ChemiDoc MP Imaging Systems Quick Start Guide, available on the Bio-Rad website.

Interacting with the Touch Screen

In Image Lab Touch, you access imaging settings by tapping the setting type on the touch screen.

- Tap an imaging settings tile to open dialog boxes in which you can specify imaging settings.
- Tap an icon to execute a command or to open a menu or a view.
- Tap a box to display an on-screen keyboard or numeric keypad in which you can enter data.

Action	Definition
Tap	Briefly touch the screen surface.
Double-tap	Tap twice quickly.
Pan	Touch and then move your finger left or right.
Swipe	Slide a finger up, down, or across a box or screen to scroll contents.
Stretch	Place a thumb and one or two fingers together on the screen, and then move them apart to zoom in.
Pinch	Place a thumb and one or two fingers slightly apart on the screen, and then move them together to zoom out.

Alternatively, you can use a USB mouse to click tiles and icons or to click in boxes, and you can enter the data with a USB keyboard.

Important: Plug in the USB mouse before you turn on the system to enable the mouse pointer to appear.

Tip: When you use a USB mouse, read the instruction *tap* in this guide as *click*.

Entering Text on the Touch Screen

You can enter or edit text in text boxes using the on-screen keyboard or keypad. You can also enter text using a USB keyboard.

To access the on-screen keyboard or keypad

- ▶ Tap in a text box.

The appropriate on-screen input device appears.

Using the On-Screen Keyboard


Text boxes appear on the screen in which you can enter or edit text.

- ▶ Tap in a box to display the on-screen keyboard.

Use the on-screen keyboard or attach a USB keyboard to the imager to enter data.

You can switch among three keyboards. Tap the keys on the bottom left of the keyboard to display the different keyboards and enter English alphabetic characters, numbers and symbols, or Western European characters.

- ▶ Choose the keyboard you want and tap keys to enter data.

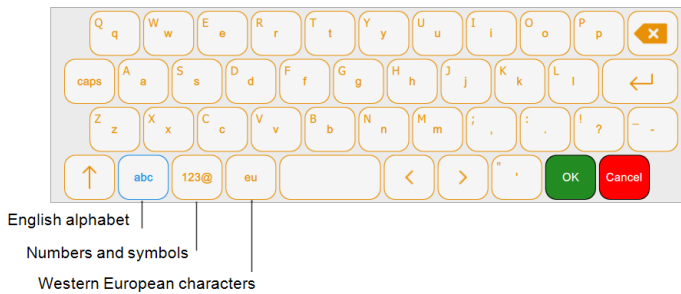


Using the On-Screen Keypad

Text boxes appear on the screen in which you can enter or edit numerals or text. You can enter numerals using the on-screen numeric keypad.

To display the numeric keypad

- ▶ Tap in a box.

The keypad appears on the screen.

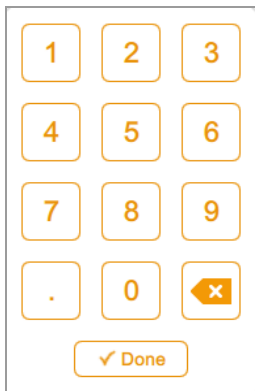


To use the keypad

- Tap the text box to display the range of valid values below the box.

- If a number you tap does not appear on the screen, verify that the number is within the range of valid values. Invalid numbers are not accepted. If you enter an invalid number and tap Done, the box turns red.
- Some properties are set by entering data in multiple text boxes. After you enter data in a box, tap the next box. All edits in each box are saved automatically.
- Tap Done or tap anywhere else on the screen to close the keypad.

Choosing Imaging Settings

Image Lab Touch displays an imaging settings tile for each channel image you want to acquire. For example, when you tap Single, Camera view displays a single imaging settings tile in the left pane:

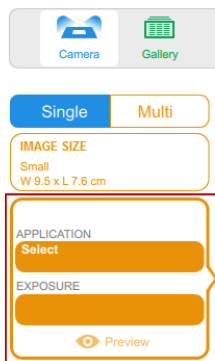

- Tap Single or Multi to acquire a single-channel or multichannel image.



- Tap Image Size to display image size settings and then select the image size you want.

IMAGE SIZE

Large

W 18.0 x L 14.4 cm

The dimensions of the image size you select appear under Image Size. Similarly, the application and exposure settings you select appear under Application and Exposure.

Close dialog box

Choose Image Size

Selected image size

Small
W 9.5 x L 7.6 cm

Medium
W 15.5 x L 12.4 cm

Large
W 18.0 x L 14.4 cm

W 18.0 L 14.4 cm

Tap to select a standard size

Tap to specify a custom size

For example, when you tap Application in the tile, application categories and application types compatible with the selected application category appear:

APPLICATION
Select

EXPOSURE
Preview

Choose Application Category

Nucleic Acid Gels

Protein Gels

Blots

Choose Application

Chemiluminescence

Colorimetric

Stain Free Blot

StarBright B700

Alexa 488

Alexa 546

Alexa 647

Alexa 680

Alexa 790

Amido Black

Cy2

Cy3

Cy5

Cy5.5

CY7

DyLight 488

DyLight 550

DyLight 650

Locking the System

If your imaging system has an administrator and if you log in with a password, you can lock the system. This prevents other users from interrupting an acquisition or changing settings.

While the system is locked, only you (the current user) or an administrator can log you off the locked system. If an administrator chooses to log you off while an action is in progress, the action and the lock are canceled.


To lock the system during an acquisition

1. Tap Camera to start an acquisition.
2. In the progress bar that appears, tap the Lock icon in the lower right corner.


The system locks and the acquisition continues.

To lock the system at any time

- On the main toolbar, tap User Settings and then tap Lock Display.

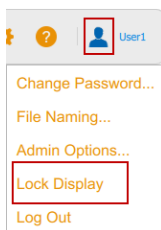

The touch screen appears dimmed. The name of the current user and buttons for unlocking or logging off the system appear in the upper right area.

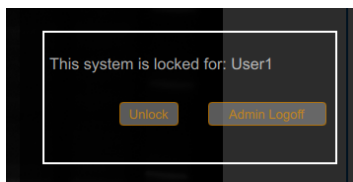

Unlocking the System

The user who locked the system (that is, the current user) can unlock it. An administrator can also unlock the system by logging off the user who created the lock. When the user is logged off, the lock is canceled, and the system is again available.

To unlock the system (current user only)


1. On the locked screen, tap Unlock.
2. Enter your password in the box that appears.

The lock is canceled and you can continue to work with the system.

To log off the current user (administrators only)

1. On the locked screen, tap Admin Logoff.

The system prompts you to identify yourself as an administrator:

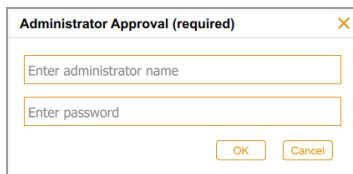A screenshot of a dialog box titled "Administrator Approval (required)" with a close button (X) in the top right corner. The dialog contains two text input fields: the first is labeled "Enter administrator name" and the second is labeled "Enter password". At the bottom right of the dialog are two buttons: "OK" and "Cancel".

Administrator Approval (required) X

Enter administrator name

Enter password

OK Cancel

2. Enter your user name and password and tap OK.

If an action is in progress, a message notifies you that data might be lost if you log off the current user.

3. Tap Yes to log off the current user or No to cancel.

Getting Help

Online help is available in each main view and in selected dialog boxes.

To access help

- Tap the question mark icon in the main toolbar to open the help topic for the current screen.
- Scroll to display longer topics and to access links to related topics.
- Tap Help Topics at the bottom of any help topic to view a list of all main help topics.

To close Help

- ▶ Tap the X in the upper right corner of the topic.

Chapter 3 Logging In to Image Lab Touch

The first time you run Image Lab Touch software, you must enter a user name to log in. You will also be prompted to indicate whether you want to be an administrator.

Logging in as a User

The first time you run Image Lab Touch software, you must enter a user name to log in. You will also be prompted to indicate whether or not you want to be an administrator.

Image Lab Touch software can be run without an administrator. In this case, users can run the software without a password and save, export, or delete their own image files. All users can update Image Lab Touch software.

To log in as a user

1. Tap anywhere in the Welcome screen.

The login dialog box appears.


A login dialog box with a light gray border. It contains a text input field with the placeholder text "Enter user name" in a light blue font. To the right of the input field is a blue plus sign icon, and below the plus sign is the word "Create" in a small blue font.

2. Tap inside the text box.
3. Using the on-screen keyboard that appears, tap in a user name and then tap Create.

The system prompts you to choose whether you want to be an administrator and displays information that helps you decide.

4. Tap No.

A Camera view of the imaging stage appears. The orange exclamation point indicates that you must take an action before you can acquire an image. Tap the exclamation point to find out more.

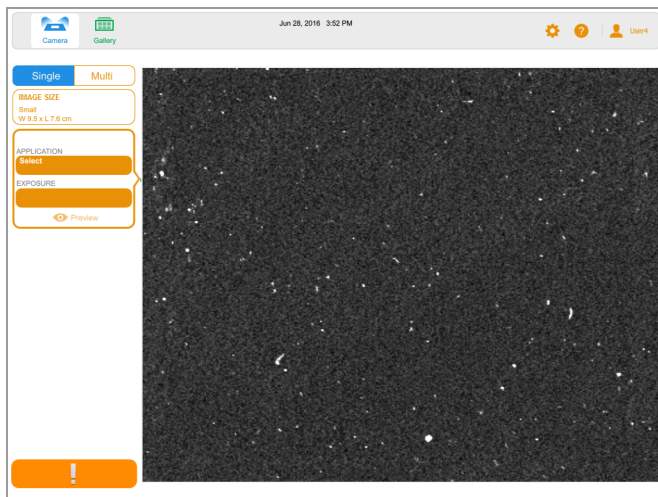

Creating a Password


After an administrator has been appointed, the administrator can require users to log in with a password and set other options to ensure the security of the system. Administrators always log in with a password.

Passwords must be 6–32 characters long and contain the following:

- one uppercase letter
- one lowercase letter
- one number
- one special character


To create a password

1. Tap User Settings in the main toolbar.

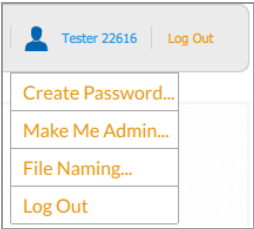

2. Tap Create Password.

The Create password dialog box appears.

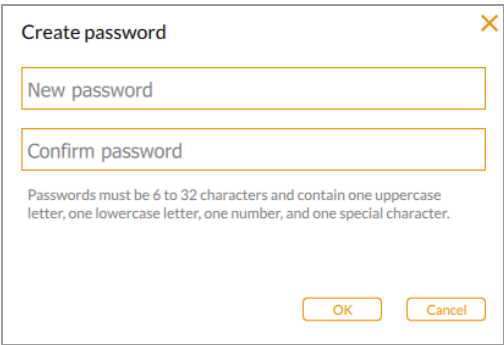

3. Tap New password.

The on-screen keyboard appears.


4. Type a password.

Tip: Tap the 123@ key in the bottom row of the keyboard to enter numbers and special characters.

5. In the Confirm password box, enter the password again.
6. Tap OK.

To change a password

1. Tap User Settings in the main toolbar.

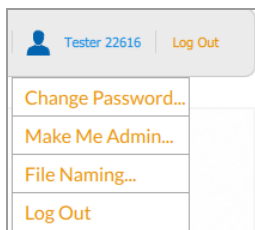

2. Tap Change Password.

The Change password dialog box appears.

A screenshot of a "Change password" dialog box. The dialog has a title bar with a close button (X). It contains three text input fields labeled "Current password", "New password", and "Confirm password". Below the fields is a note: "Passwords must be 6 to 32 characters and contain one uppercase letter, one lowercase letter, one number, and one special character." At the bottom right are "OK" and "Cancel" buttons.

3. Tap Current password.

The on-screen keyboard appears.

4. Type your current password.
5. Enter a new password in the New password box.
6. Enter the new password again in the Confirm password box.
7. Tap OK.

Removing a Password

If no administrator has been assigned or if the administrator does not require users to log in with a password, you can remove your password.

To remove a password

1. Tap User Settings in the main toolbar.
2. Tap Change Password in the dropdown menu.
3. In the Change Password dialog box, delete your current password and tap OK.

Logging in as an Administrator

Any user can become an administrator at any time. However, once an administrator has been appointed, only the administrator can create other administrators.

Tip: Store your administrative password somewhere safe. For security purposes, administrative password recovery requires a field service visit.

To log in as an administrator


1. Tap anywhere in the Welcome screen.
2. Using the on-screen keyboard that appears, tap in a user name and then tap Create.

The system prompts you to choose whether you want to be an administrator and displays information that helps you decide.

3. To become an administrator, tap Yes and create a password in the box that appears.

The Administrator Options dialog box appears, in which you can choose options for managing Image Lab Touch user accounts.

Chapter 3 Logging In to Image Lab Touch

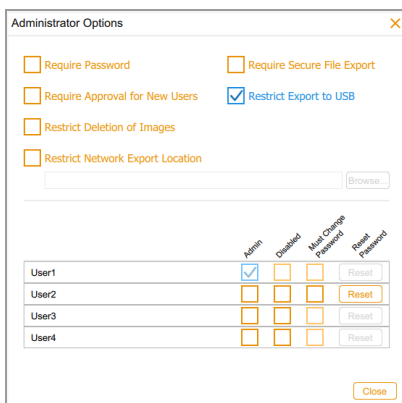

The Administrator Options dialog box contains several settings and a user management table.

Settings:

- ☐ Require Password
- ☐ Require Secure File Export
- ☐ Require Approval for New Users
- ☒ Restrict Export to USB
- ☐ Restrict Deletion of Images
- ☐ Restrict Network Export Location

User Management Table:


	Admin	Disabled	Max Change Password	Reset Password
User1	<input checked="" type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>
User2	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>
User3	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>
User4	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>

Buttons: Close

See [Table 1 on page 46](#) for information about the options in this dialog box.


4. Choose the options you want and tap Close.

A Camera view of the imaging stage appears.

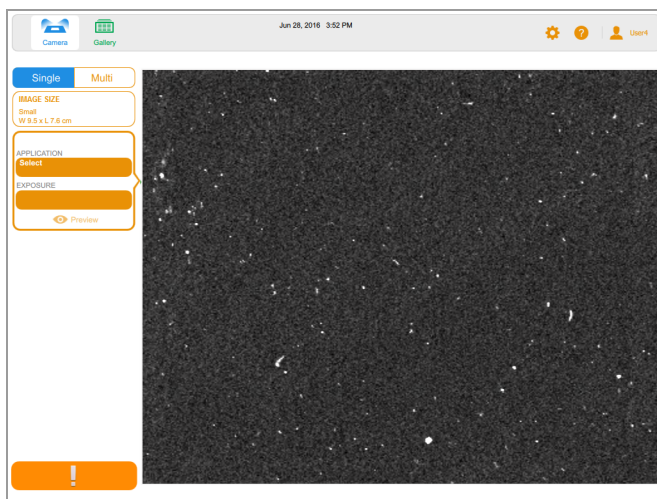

To become an administrator after login

1. In any view, tap Make Me Admin in the User menu.

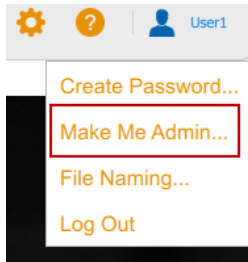

The Administrator Options dialog box appears, in which you can choose options for managing Image Lab Touch user accounts.

Setting Administrator Options

Having an administrator adds security and file access control features to the imaging system. When you log in as an administrator, you can set the following administrator options.

Table 1. For All User Accounts

This Option	When Selected
Require Password	Requires all users to log in with a password. If the user has no password, they must create one.
Require Approval for New Users	Requires administrator to approve all new user accounts.
Restrict Deletion of Images	Prevents all users who are not administrators from deleting images.
Require Signed File Export	<p>Requires files to be exported as secured documents that cannot be altered without detection and that can be used as protected documents in Image Lab, Security Edition.</p> <p>Note: Selecting this option also enables the options to require passwords and to make new user accounts subject to administrator approval.</p>
Prohibit Export to USB	Prevents users from exporting files via USB.
Restrict Network Export Location	Restricts users to a top-level network location to which they may export files.

Table 2. For Individual User Accounts

This Option	When Selected
Admin	Makes this user an administrator.
Disabled	Disables this user's account. Note: Only administrators can access disabled accounts.
Must Change Password	Requires this user to reset password.
Reset Password	Allows the administrator to enter a new password for a user.

To set administrator options

1. Tap User Settings and then tap Admin Options.

The Administrator Options dialog box appears.

Administrator Options

☐ Require Password

☐ Require Secure File Export

☐ Require Approval for New Users

☒ Restrict Export to USB

☐ Restrict Deletion of Images

☐ Restrict Network Export Location

	Admin	Disabled	Must Change Password	Reset Password
User1	<input checked="" type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>
User2	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>
User3	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>
User4	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="button" value="Reset"/>

2. Select the checkbox for each option you want to enforce.
3. Clear the checkbox for each option you want to allow.

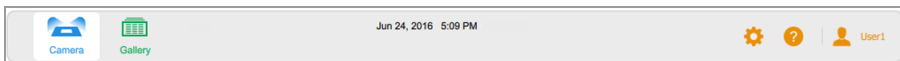
Chapter 4 Understanding the Views

Image Lab Touch software guides you through the processes of acquiring and managing images in four views. Color Cues in these views signal different states in the system.







- Camera View
- Preview
- Image View
- Gallery

Main Toolbar Settings

The main toolbar appears at the top of every Image Lab Touch view.




Tapping a toolbar object opens a view or a menu.

Object	Explanation
	Camera view — displays a view of the sample in the sample tray in real time. Also displays a preview image of the sample.
	Gallery — displays thumbnails of all images you acquired. You can view, browse, delete, print, or export images. Double-tapping a thumbnail displays the image full size in Image View.
	Retake — visible when you acquire a multichannel image, enables you to retake one or more channel images right after acquisition.
[Date and time]	Date and Time — displays the current date and time. To change this setting, go to System Settings > Set Date and Time.
	System Settings — menu that displays system settings and accesses software updates. Also accesses white tray calibration settings when a white tray is inserted for the first time.
	Help — displays procedures and concepts for the view on screen.
	User Settings — menu that accesses commands to change your password, set a file naming format for acquired images, view Admin settings, lock the display, or log out. Name of current user appears to its right.

Camera View


When you log in to the imaging system, Camera view displays the imaging stage.

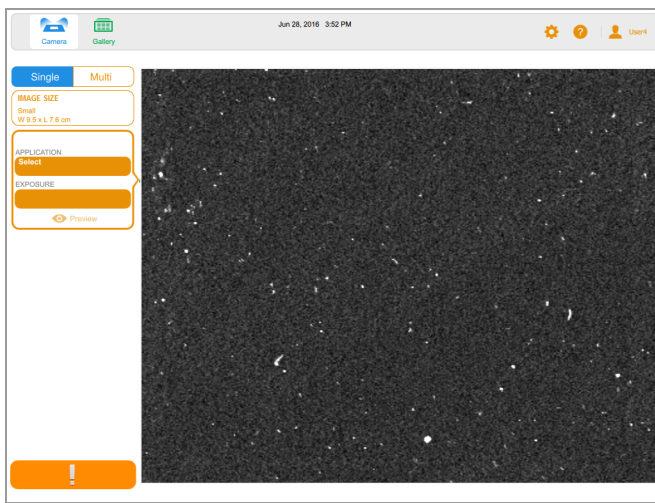

When you place a sample in the imager, Camera view displays the sample in real time. The left pane displays imaging settings tiles for each channel you want to acquire. You can specify an image size, select an application, and set the exposure for each channel.

After you choose these settings, you can acquire the image.

The acquired image appears full size in Image View and also as a thumbnail in the Gallery.

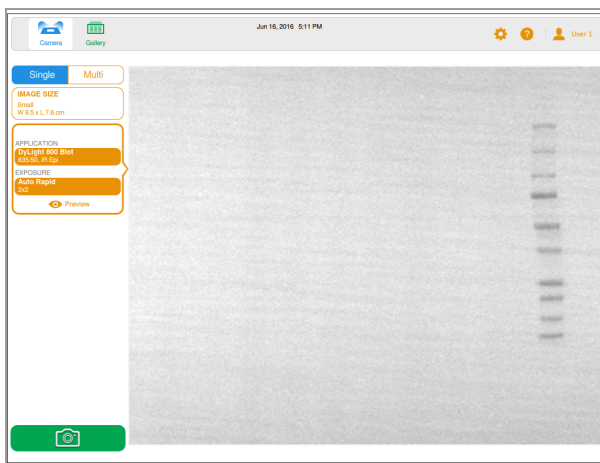

- **Main toolbar** — accesses Camera view, the Gallery, help topics, system settings, and user settings. For more information, see [Main Toolbar Settings on page 49](#).
- **Imaging area** — displays the sample in the imager.
- **Image Settings Tiles** — tapping Application or Exposure in a tile displays options for that setting.
- **Preview** — tapping Preview in the tile displays a low-resolution image of the sample, which you can use to assess the image.
- **Camera button** — tapping this button at the bottom left of the screen acquires an image of the sample.

For more information, see [Acquiring an Image — General Steps on page 74](#).

Preview

Preview quickly displays a low-resolution image of the sample in the imager so you can determine whether sufficient sample is visible and identify the region for which you want to set an automatic exposure.



For more information, see [Previewing an Image on page 81](#).

Image View

When you acquire an image, Image View displays it full size. You can also select up to four image thumbnails at a time in the Gallery. The selected images open full size in Image View.



Settings in the Image View toolbar at the bottom of the screen enable you to fine tune and manage images.

The Image View Toolbar


The Image View toolbar appears at the bottom of Image View.



Icon	Description
	Visible when a single image is displayed. Tap to scroll through Previous and Next images selected for display in the Gallery.
	Copy Zoom — visible when multiple images are displayed. Selecting a zoomed image and then tapping this icon copies the zoom setting of the selected image to the other displayed images.
	Transform — enables you to optimize image appearance by adjusting brightness and contrast.
	Image Info — lists acquisition details for the active image and the name of the user who acquired it. You can rename an image or add notes about it.
	Merge — combines up to three images into a multichannel image.
	Print — prints the displayed image to the local printer.
	Delete — deletes the displayed image.
	Send/Save — saves the displayed image to a USB flash drive or network drive.

The Gallery


The Gallery displays thumbnail images of all acquired images in order by date acquired, most recent acquisition date first. Each thumbnail image is named with the date and time of acquisition. You can change this name in the Image Info box.

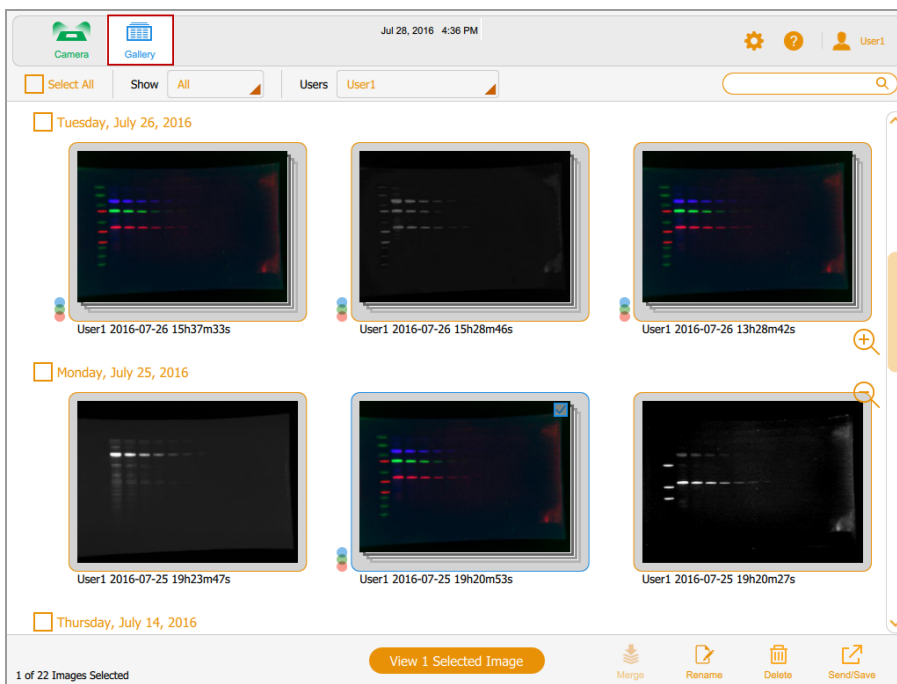

Tap a thumbnail image in the Gallery to open it full size in Image View, delete selected images, and save selected images to a USB flash drive, an external hard drive, a stand-alone computer, or a network drive. Tap zoom icons to make thumbnails bigger or smaller.

For more information, see [Chapter 8, Viewing Images](#).

To compare images, you can select up to four images in the Gallery to display in Image View.

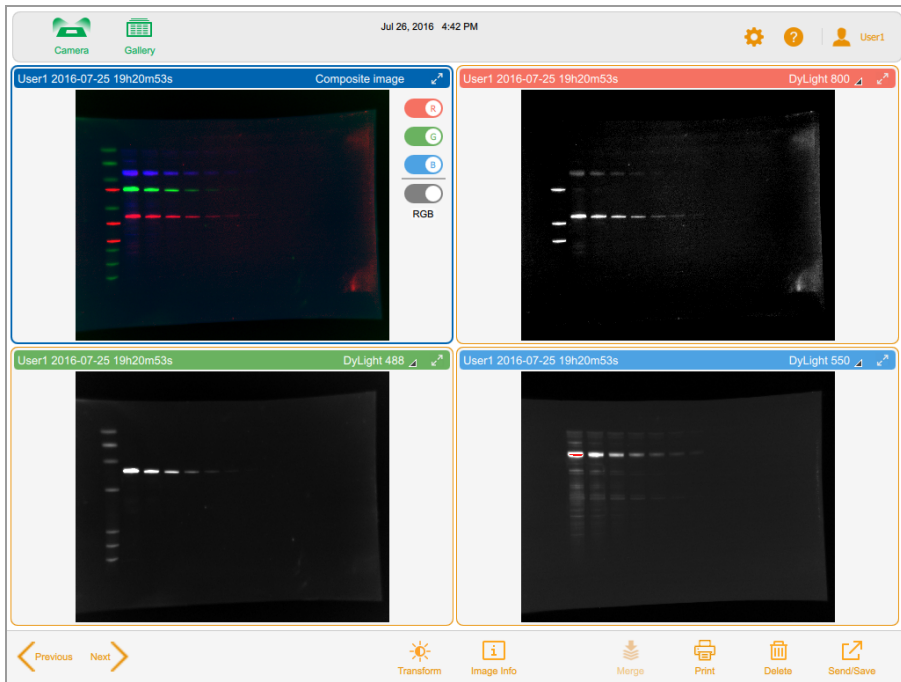


For more information, see [Comparing Images on page 123](#).

Color Cues


Image Lab Touch color cues signal you to the various states of the system.

Note: In this section, the definition of each color cue is followed by an example screen, in which a red outline highlights the use of the color cue.

Green and **Orange** indicate active objects. Here, the green color cue indicates that you can select Gallery. The orange exclamation point indicates that an error message requires action.

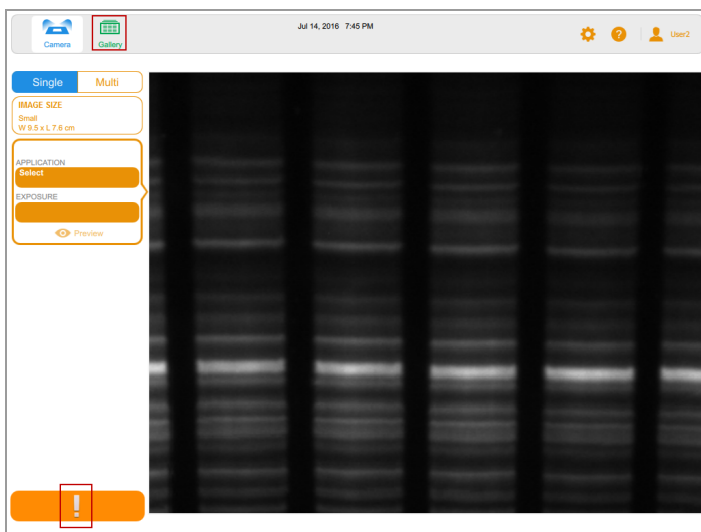


Blue indicates a selected object. In this image, Camera view and single image are selected.

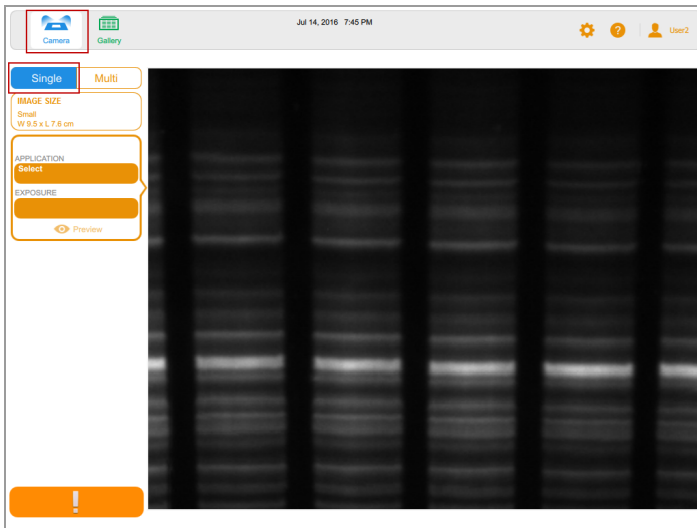


In the Gallery and in Image View, selected images are outlined in blue.


Chapter 4 Understanding the Views

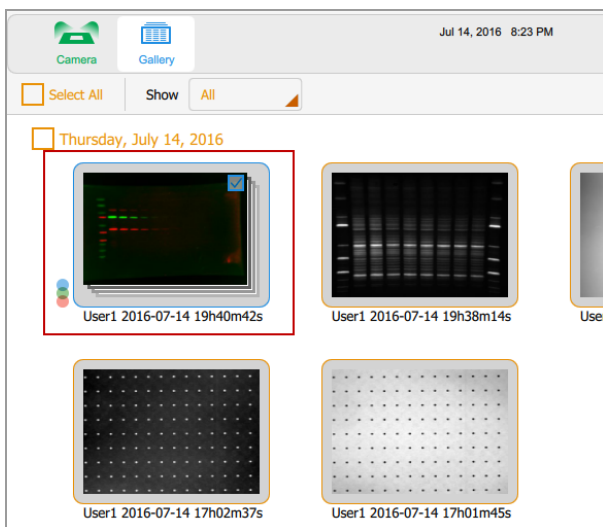


Orange indicates that an action is required before you can continue. In the next image, the orange exclamation point appears in place of Camera to indicate that an error message requires action.

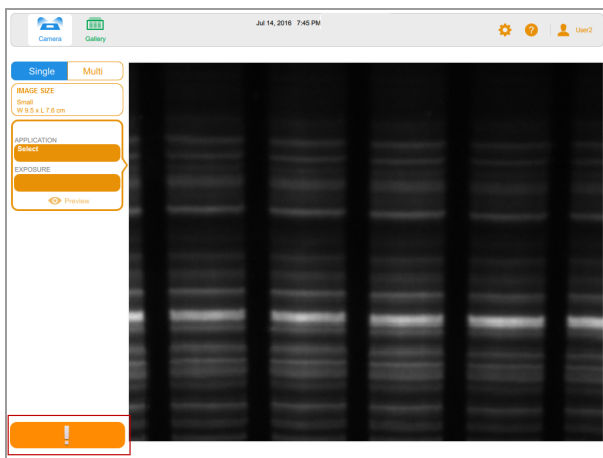

Tap the orange exclamation point to display a message that explains what to do.

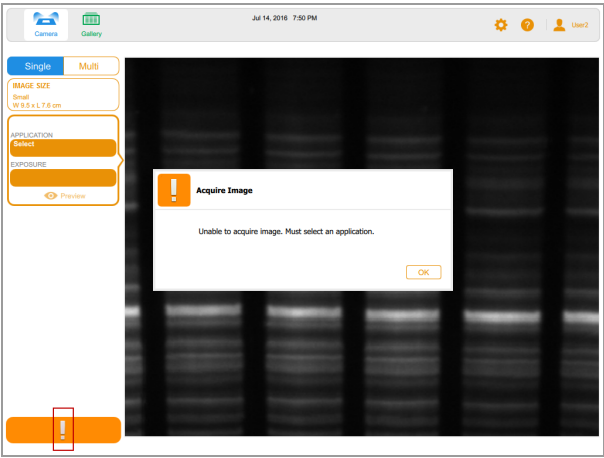

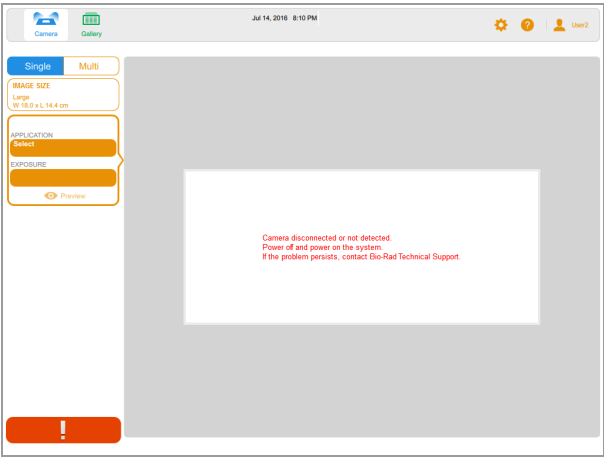


Red indicates a critical error. Tap the red exclamation point to display the message. You must log out of Image Lab Touch and restart the imager to correct the error.



Chapter 5 Choosing a Sample Tray

Before you acquire an image, select a sample tray designed for the application you select.


The imaging system supports the following sample trays:

Sample Tray Name	Icon
Blot/UV/Stain-Free tray	
White tray	
Blue tray	

The imager detects the type of sample tray on the imaging stage. If the tray in the transilluminator drawer does not support the application, an error message appears and the imager does not acquire the image.

Some fluorescent reagents have excitation peaks in both UV and blue wavelengths. In this case, either excitation source can be used. However, the resulting images are not identical. Using the recommended tray with these fluorescent reagents provides a better signal-to-noise ratio, which increases sensitivity.

Tip: When two tray types can be used, both are shown. An asterisk identifies the recommended tray.

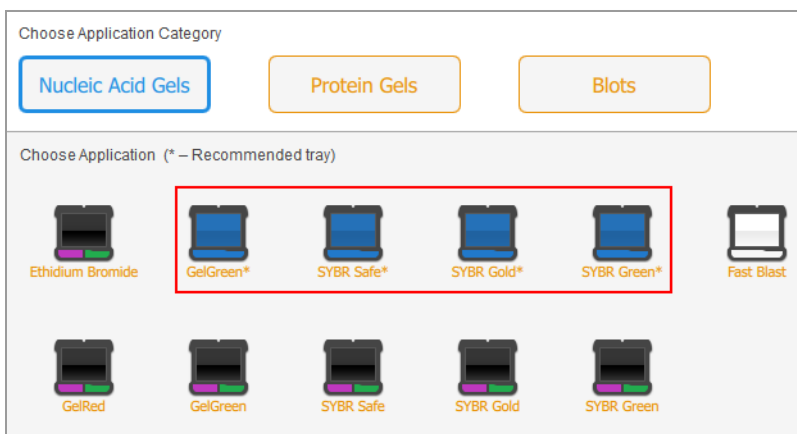

To position a sample on the sample tray

1. Place the sample face up and center it on the appropriate sample tray.

Tip: Center the sample carefully so it fills the image area. If you enlarge the image by zooming in, check to ensure that the sample is still centered.

2. Open the imager door and pull out the transilluminator drawer.
3. Pull out the transilluminator drawer.
4. Place the sample tray on the imaging stage.
5. Close the transilluminator drawer and the door.

[Applications Listed by Tray Type](#)

Applications Listed by Tray Type














The ChemiDoc and ChemiDoc MP Imaging Systems ship with the Blot/UV/Stain-Free Tray. Two optional trays are available for purchase. Applications you can use with each tray are listed in the following tables.

Blot/UV/Stain-Free Tray

Some applications for this tray can be processed on both ChemiDoc and ChemiDoc MP Imaging Systems for certain applications. Others can be processed only on ChemiDoc MP, as indicated in the following table.

Applications for the Blot/UV/Stain-FreeTray (cat. #12003028)		
	ChemiDoc	ChemiDoc MP
Ethidium Bromide	■	■
GelStar	■	■
Krypton	■	■
Ponceau S	■	■
Pro-Q Diamond	■	■
Pro-Q Emerald 300	■	■
Pro-Q Emerald 488	■	■
GelGreen	■	■
Stain-free gel	■	■
Stain-free blot	■	■

Applications for the Blot/UV/Stain-FreeTray (cat. #12003028)		
	ChemiDoc	ChemiDoc MP
Colorimetric	■	■
Chemiluminescence	■	■
Amido Black	■	■
SYBR® Gold	■	■
SYBR® Green	■	■
SYBR® Safe	■	■
StarBright B520*	■	■
GelRed	■	■
Rhodamine		■
StarBright B700		■
IRDye 800		■
IRDye 680		■
Alexa 488		■
DyLight 800		■
Alexa 546		■
Alexa 647		■
Alexa 680		■

Applications for the Blot/UV/Stain-FreeTray (cat. #12003028)		
	ChemiDoc	ChemiDoc MP
Alexa790		
Coomassie Blue**		
CY2		
SYPRO Ruby blot		
CY3		
CY5.5		
CY7		
DyLight 488		
DyLight 550		
DyLight 650		
DyLight 680		
CY5		
SYTO 60		

* For best results image StarBright B520 on ChemiDoc MP.

** For best results image Coomassie Blue using the blot/UV/stain-free tray. Doing so allows the Coomassie Blue dye to excite with Far Red light, which produces a signal to noise ratio superior to traditional white light.

Blue Tray

Applications for the Optional Blue Tray (cat. 12003026)

- | | |
|--------------|---------------|
| ■ GelGreen | ■ SYBR® Green |
| ■ GelStar | ■ SYBR® Safe |
| ■ SYBR® Gold | |

White Tray

The white tray must be calibrated before it is used for the first time.

Applications for the Optional White Tray (cat. 12003026)

- | | |
|------------------|----------------|
| ■ Coomassie Blue | ■ Silver stain |
| ■ Fast Blast | |

Calibrating the White Tray


The optional white tray requires calibration before it can be used for the first time.

Note: If you subsequently obtain another white tray, Bio-Rad recommends that you calibrate the new tray.

When the imaging system detects that a white tray has been inserted for the first time, a message indicates that the tray is not calibrated and lists steps to calibrate it.


To calibrate the white tray

1. Complete the steps listed in the message box.

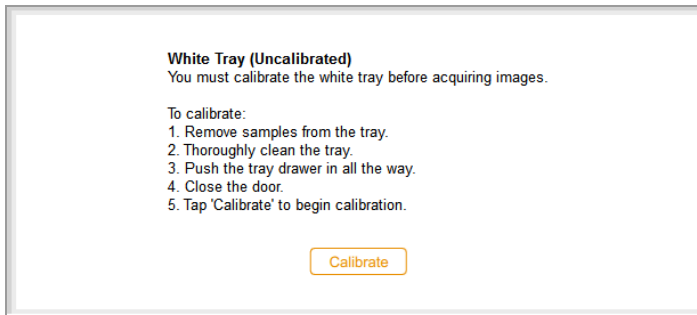

2. Tap Calibrate.

The system starts calibrating the white tray. A progress bar monitors the process.

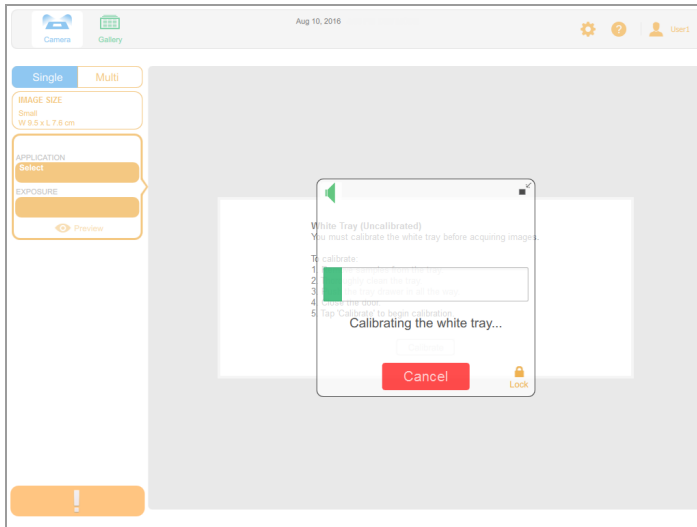

A message notifies you when calibration is complete.


3. Tap OK to close the message.

Calibrating a Replacement White Tray


When you replace the white tray, Bio-Rad recommends that you calibrate the new tray.

To calibrate a replacement white tray

1. Tap User Settings, and then tap Calibrate White Tray.

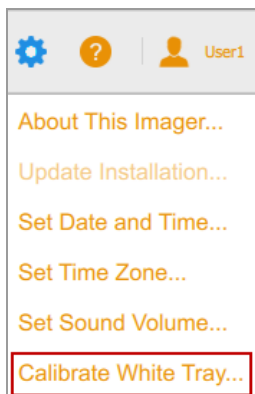

A message lists steps to calibrate the tray.

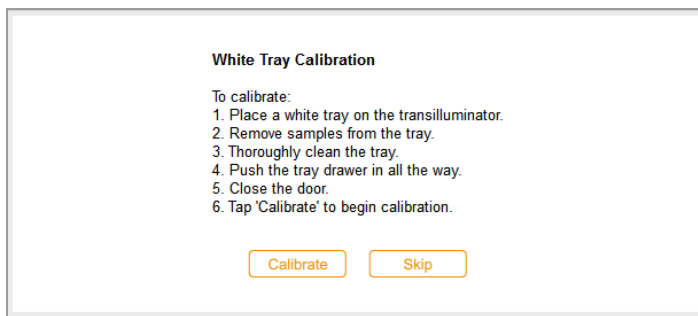

2. Complete the steps listed in the message box.
3. Tap Calibrate.

The system starts calibrating the white tray. A progress bar monitors the process.

A message notifies you when calibration is complete.

4. Tap OK to close the message.

Chapter 6 Acquiring an Image

Instructions for acquiring an image apply to both single- and multichannel images unless they are preceded by (Single Channel) or (Multichannel).

The steps for acquiring single-channel and multichannel images are similar. Acquiring a multichannel image requires repeating some steps for additional channels, and a few procedures are specific to single- or multichannel images.

Before you can acquire an image, you must place a sample on the sample tray and then place the tray on the imaging stage. You can specify acquisition settings before you place the sample on the sample tray or after you place the tray in the transilluminator drawer. The imager retains your settings until you change them.

Multichannel Images

Note: Only applications for blots are available for multichannel images.

You can configure 1–3 channels on the ChemiDoc MP and 1–2 channels on the ChemiDoc imaging system for multichannel acquisition. The image size you choose applies to all channels in the acquisition.

When you choose an application for a channel, a plus sign surrounded by a dotted outline tile appears below the selected channel. Tapping inside the outline box changes it into an imaging settings tile for the next channel.



Placing the Sample in the Imager

You must use the tray that supports the application you select. The imaging system detects the type of sample tray on the imaging stage. If the tray in the transilluminator drawer does not support the selected application, an error message appears and you will not be able to acquire the image.

To place the sample in the imager

1. Place the sample face up on the sample tray.
Tip: If you plan to zoom in or out on the image, center the sample on the tray as accurately as possible.
2. Open the imager door.
3. Pull out the transilluminator drawer.
4. Place the sample tray on the imaging stage.
5. Close the transilluminator drawer and the image door.

Acquiring an Image — General Steps


General steps outline the workflow for acquiring an image. Step-by-step procedures in the following sections guide you through these general steps.

Note: You can configure 1–3 channels on the ChemiDoc MP and 1–2 channels on the ChemiDoc Imaging System for multichannel acquisition. The image size you choose applies to all channels in the acquisition.


To acquire an image

1. In Camera view, do one of the following:

- (Single Channel) Tap Single to configure a single-channel image.

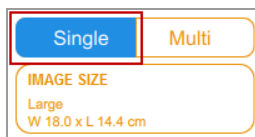


- (Multichannel) Tap Multi to configure a multichannel image.

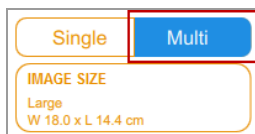

2. Specify the image size so that the sample fills as much of the screen as possible.
3. (Multichannel) Select a channel color or keep the default color. (You can change the channel color at any time.)
4. Choose the application.
5. Preview the image.
6. Keep the default automatic exposure option or change the exposure options.
Note: When you set an automatic exposure time and specify a region of interest, the optimal auto-exposure time applies to that region only.
7. (Multichannel) Repeat steps 3–6 for the remaining channels.
8. Acquire the image.

Setting Image Type and Size

Image type and size settings appear at the top of the left pane in Camera view. In multichannel acquisitions, the image type and size you choose apply to all channels.

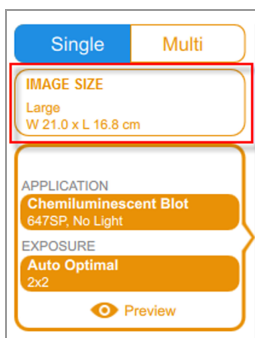

Choosing Image Type

To choose an image type

- ▶ Tap Single or Multi to acquire a single-channel or multichannel image.



Choosing Image Size


You can choose a preset image size or specify a custom size. The preset sizes correspond to these Bio-Rad gels:

- **Small** — Mini-PROTEAN
- **Medium** — Criterion
- **Large** — Wide Mini ReadyAgarose

To choose a preset image size

- ▶ Tap Image Size to select a size in the dialog box that appears.

The dimensions of the selected image size appear under Image Size.

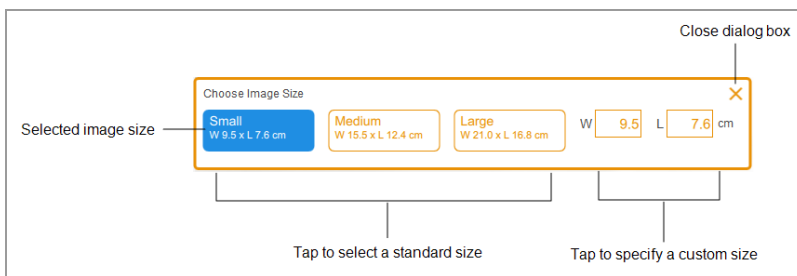

To specify a custom image size for a gel or blot image

1. In Camera view, tap Image Size.
2. Tap the W (width) or L (length) box. When you enter one dimension, the software automatically calculates the other according to the imager's 5:4 aspect ratio.
3. Using the keypad, enter the width or length of the image (in cm).

Valid width values are 9.0–21.0 cm. Valid length values are 7.2–16.8 cm.

To change the image size manually

- ▶ Place a thumb and one or two fingers slightly apart on the screen and move them together (pinch to zoom) without lifting them from screen until the image is the size you want.


Choosing the Application

You can choose from a list of predefined applications for common sample types.


Image Lab Touch software supports three application categories: nucleic acid gels, protein gels, and blots. Multichannel images can be acquired only for blots.

Important: You must use the tray that supports the application you select. The imager detects the type of sample tray on the imaging stage. If the tray in the transilluminator drawer does not support the selected application, an error message appears when you try to acquire the image.

When you choose an application, the software selects the appropriate filter and illumination source. Detailed settings appear in the imaging settings tile under the name of the selected application. The default automatic exposure setting also appears. You can change the default exposure setting.

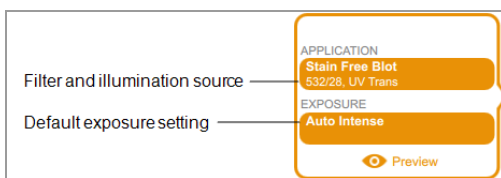


When you select a low-light application, the optimal binning level also appears under Exposure.

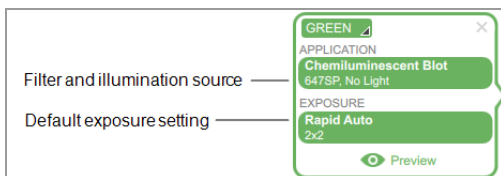

To choose an application

1. In Camera view, tap Application in an imaging settings tile.


The Application box appears.

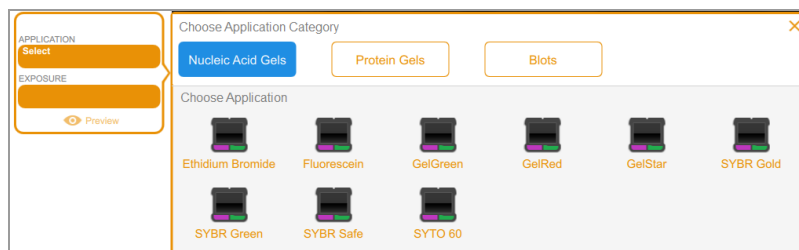

2. Under Choose Application Category, tap an application category.

The Choose Application list displays only applications compatible with the category you selected.

For example, when you tap Blots, only applications for blots appear.

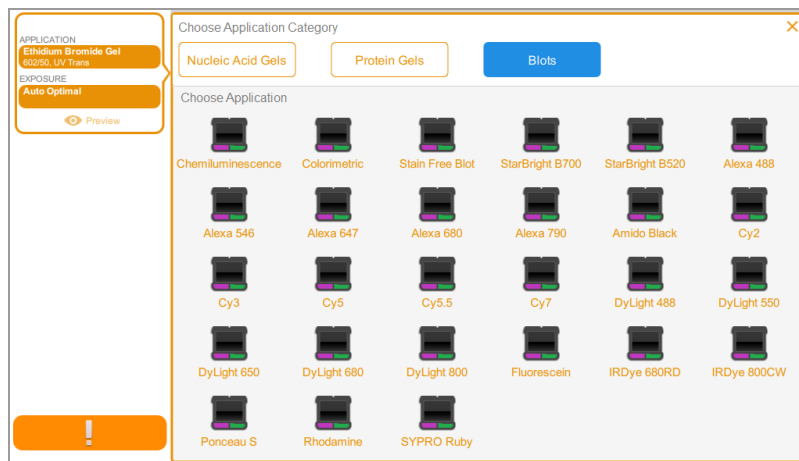

3. Tap an application.

(Multichannel) When you configure the first channel, the system eliminates applications that are redundant or incompatible with the application you selected, and so on for the second and third channels you configure.

4. (Multichannel) Repeat steps 1–3 for each channel you want to add.

(Multichannel) Selecting a Channel Color


Assigning channels different colors makes it easier to identify the channels in a composite image. You can use any channel color with any application when you specify channel settings.

Note: There is no relationship between the emission wavelength of the sample and the channel color.


Acquisition options for a channel display a default channel color. You can keep the default color or change it. You can change a channel color at any time. If you change the channel color to one already assigned to another channel, the channels swap colors.

To select a channel color

1. Tap to view the color list.

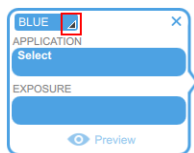


2. Tap a color to select it.

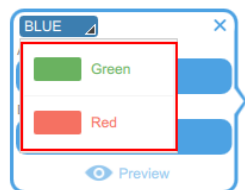

Previewing an Image

When you put a sample in the imager and close the door, Preview appears at the bottom of the imaging settings tile.

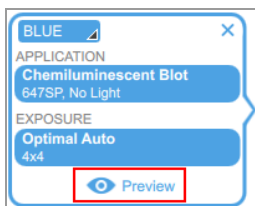


When you tap Preview, the system displays a quick, low-resolution image of the sample using the appropriate light source and filter for the application. View this image to determine whether sufficient sample is visible. For an auto exposure, you can also identify a region of interest. If sufficient sample is visible, you can then select either Auto Optimal or Auto Rapid exposure for the image you want to acquire.

Important: Keep the imager door closed while viewing the preview image.

To preview an image

1. Specify an application.
2. Tap Preview to view a low-resolution image of the sample.

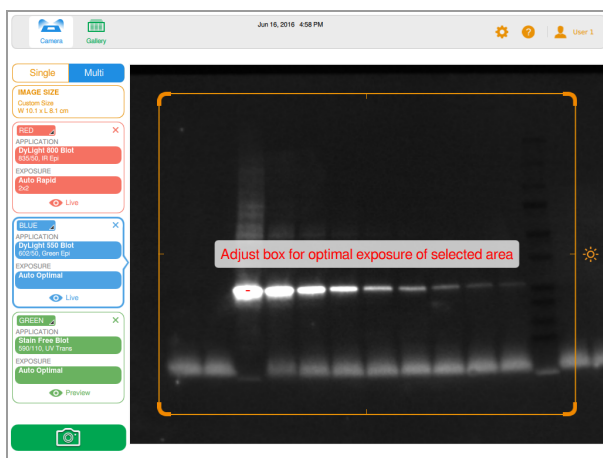

3. Inspect the preview image to verify that important sample features are visible in the sample.
4. (Optional) For an automatic exposure, specify a region of interest (for example, a lane or band).
5. Tap Camera to start the Auto Exposure for the defined region of interest.

To clear the preview image, do any of the following:

- Open the door.
- Select a different application.
- Switch between single and multichannel modes.
- Log out.
- Delete a channel and then add it again.

Exposure Settings

Steps for specifying exposure settings are specific to the type of application you select and how you choose to expose it: automatically (Rapid or Auto) or manually.

Refer to the section about the type of application you want to image.

- Most gels and blots
- Stain-free gels
- Chemiluminescent and low-light blots

Exposure Settings for Most Gels and Blots

To choose an automatic exposure setting for a gel or blot

1. In Camera view, tap Exposure.
2. Tap Auto.
3. Tap Optimal or Rapid Auto-exposure.
4. (Optional) Tap Preview and then specify the region of interest in the preview image.
5. Tap Camera to acquire the image.

To manually set an exposure time for a gel or blot

1. In Camera view, tap Exposure.
2. Tap Set Manual Exposure Time.
3. Enter the time in the text box using the numeric keypad.
4. Tap Camera to acquire the image.

[Exposure Settings for Chemiluminescent and Low-Light Blots](#)

Exposure Settings for Stain-Free Gels

Before you set the exposure time for a stain-free gel, choose a gel activation time.

To set the gel activation time

1. Tap Exposure.
2. Under Select Gel Activation, choose a gel activation time:
 - **No Activation** — to retake an image for a previously activated gel.
 - **45 sec** — if you plan to transfer the gel to a blot for immunodetection.
 - **5 min** — when you want to detect proteins that are low in concentration and you plan to quantitate the maximum number of bands. This duration provides an optimal signal-to-noise ratio because gel activation is close to completion at the end of five minutes.

To choose automatic exposure settings

1. Tap Auto.
2. Choose Optimal or Rapid Auto-exposure.

Tip: When you choose Optimal, the dynamic range of the image is maximized to show faint bands without saturating the bright bands. When you specify a region of interest, the dynamic range is optimized for the specified area only.

Tip: When you choose Rapid, the software calculates a reasonable exposure time more quickly than Optimal. The resulting image might have less dynamic range, and intense bands and other features might be saturated.

3. Tap Camera to acquire the image.

To set a manual exposure time

1. Tap Manual.
2. Tap the sec (seconds) box to display the keypad.
3. Enter the exposure time in seconds.
4. Tap Done.
5. Tap Camera to acquire the image.

Setting Exposure Time for a Chemiluminescent or Low-Light Blot

You can choose automatic or manual exposure settings and adjust the binning level for a chemiluminescent or low-light fluorescence application (Far/Infra Red).

Low-light applications include the following:

- CY5.5
- CY7
- Alexa 680
- Alexa 790
- DyLight 680
- DyLight 800
- IRDye 680RD
- IRDye 800CW

Image Lab Touch software estimates the best exposure time for obtaining good-quality images with the optimal dynamic range from low-light applications.

Automatic Exposure Settings

You can choose how Image Lab Touch calculates an automatic exposure. Two automatic exposure settings are available:

- **Optimal** — for an image that maximizes the dynamic range to show faint bands without saturating the bright bands. When you specify a region of interest, the dynamic range is optimized for the selected area only.
- **Rapid** — for a quicker estimate of the exposure time than Optimal. The exposure itself might not take less time, but the software calculates a reasonable exposure time more quickly. The resulting image might have less dynamic range, and intense bands and other features might be saturated.

Tip: Using Rapid Auto-exposure can help you estimate manual exposure time. Choose Rapid Auto-exposure and acquire the image. When the acquired image appears, tap Image Info in the lower toolbar to view the exposure time. Using this exposure time as a baseline, manually set the exposure time to obtain the image quality you want.

Setting Exposure Automatically for Chemiluminescent or Low-Light Applications

Setting Exposure Automatically for Chemiluminescent or Low-Light Applications


You can choose automatic or manual exposure settings and adjust the binning level for a chemiluminescent or low-light fluorescence application (Far/Infra Red). Image Lab Touch software estimates the best exposure time for obtaining good-quality images with the optimal dynamic range from low-light applications.

By default, the software selects a binning level based on either the selected application or on a combination of application and image size.

To choose an automatic exposure setting for a chemiluminescent or low-light application

1. In Camera view, tap Exposure.
2. Review the recommended binning setting on the scale and tap a different binning setting if you like.

Tip: Bio-Rad recommends starting with the optimal binning setting, examining the result, and then adjusting settings when acquiring subsequent images.

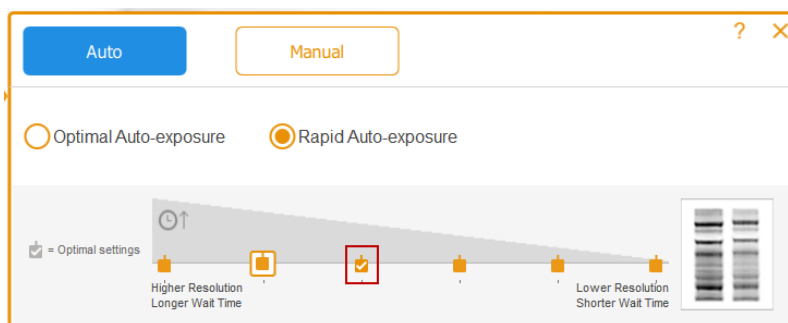

Legend

- Binning level
- Selected binning setting
- Optimal binning setting

3. Tap one of the following:
 - Optimal Auto-exposure
 - Rapid Auto-exposure

4. (Optional) Tap Preview to obtain a preview image and then specify the region of interest.
5. Tap Camera to acquire the image.

Specifying the Region of Interest for Auto-Exposure

Specifying a region of interest is especially valuable when setting automatic exposure for chemiluminescent or fluorescent samples.

Note: You cannot specify a region of interest for a stain-free gel.

When you preview an image set for auto-exposure, you can specify a region of interest such as a band. Image Lab Touch sets the automatic exposure so that the brightest pixels in the region of interest are exposed to near saturation.

Tip: Because protein standards are intense, consider excluding them from the region of interest. If you do not exclude standards, fainter features might not be visible.


Consider specifying the region of interest when you want to do the following:

- To evaluate relative signal intensities, all features that you want to compare must be unsaturated. To prevent saturation of features, specify a region that encompasses the features to be compared.
- To push the limits of sensitivity, specify a region of faint or no apparent signal. The software optimizes the exposure to reveal features in the specified region, but this might cause saturation of more intense features on the blot.

Important: Keep the door closed during preview. Opening the door clears the preview image.

To specify the region of interest in a preview image


1. Tap Preview in the imaging tile and wait for the light source to warm up.

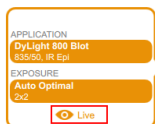

A low-resolution image of the blot appears with a selection rectangle you can adjust to surround an area of interest.




2. To resize the rectangle, drag a corner.
3. To move the rectangle, drag an edge.

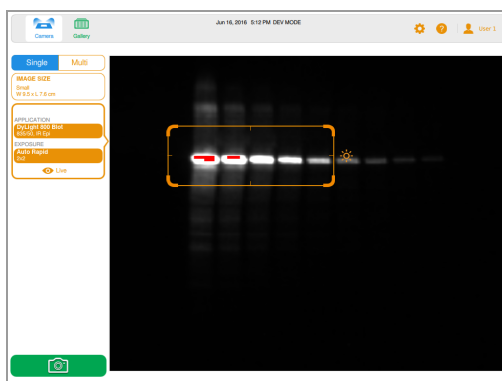

4. To acquire the image, tap Camera.

In Camera view, Preview now reads Live.

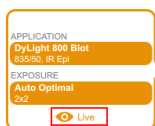

5. Tap Live to return to a view of the sample on the imaging stage.

Setting the Exposure Manually


For a manual exposure time, the following options are available:

- **Set Manual Exposure Time** — to enter time manually in seconds.
- **Configure Signal Accumulation Mode** — (chemiluminescent blots and single-channel only) to have the imager capture a series of images with a range of exposure times.




To choose a manual exposure setting for a chemiluminescent or low-light blot

1. In Camera view, tap Exposure.
2. Tap Manual.

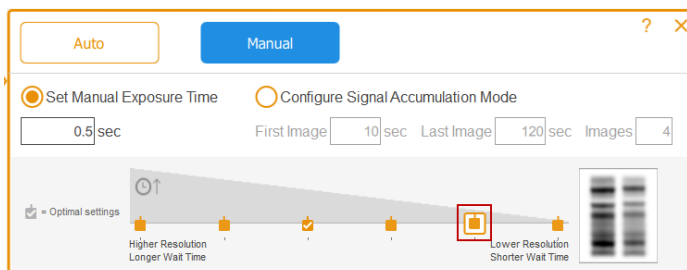- Review the recommended binning setting on the Image Resolution/Time scale and tap a different binning setting if you like.



Legend

-  Binning level
-  Selected binning setting
-  Optimal binning setting

- Tap Set Manual Exposure Time.
- Tap the sec (seconds) box to display the keypad.
- Enter the exposure time in seconds.

Tip: You can use the exposure time from a Rapid Auto-exposure image as a baseline for estimating exposure time.

- Tap Camera to acquire the image.

(Single Channel) Acquiring Signal Accumulation Mode (SAM) Images

SAM is useful for determining the best exposure time for a chemiluminescent sample.

SAM produces a series of cumulative images acquired using the range of exposure times you set. After using SAM to determine the best exposure time, Bio-Rad recommends that you acquire a single image of the sample using the optimal exposure to reduce the amount of noise added by the accumulation process.

During SAM acquisitions, a preview of each image appears so you can monitor its quality.

To calculate the SAM exposure setting

- Estimate the shortest and longest exposure times you think will achieve the optimal image.



For example, if you think the minimum time to acquire an image is 60 seconds and the longest time is 180 seconds, set these as the first and last values and then define the number of images to be acquired during the interval.

Tip: If no features appear in the resulting SAM images, acquire a second group of SAM images using longer exposure times.

To acquire SAM images

1. In Camera view, tap Exposure.
2. Tap Manual and then tap Configure Signal Accumulation Mode.
3. Enter the exposure times, in seconds, in the First Image and Last Image boxes.

(Single Channel) Acquiring Signal Accumulation Mode (SAM) Images

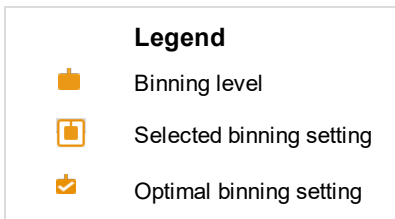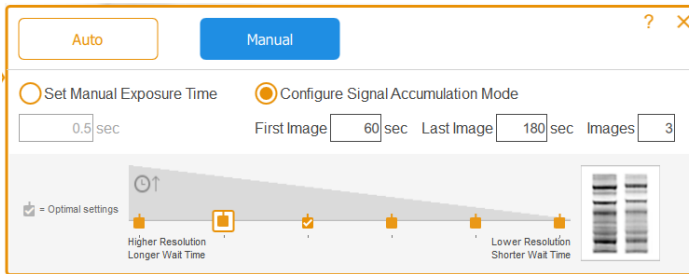


4. Enter the total number of images in the Images box. The recommended total is 3–5 images.
5. Tap Camera to acquire the image.


The last image acquired appears full size in Image View.

To obtain the exposure time for the preferred SAM image


1. In the Gallery, tap the SAM group icon.

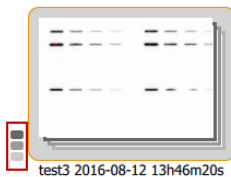

Individual SAM image thumbnails appear.

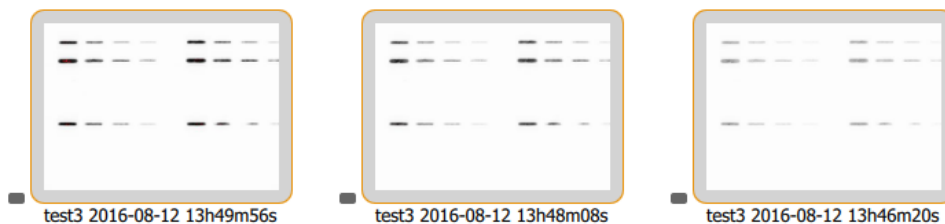

2. Tap the image with the exposure you prefer to select it.

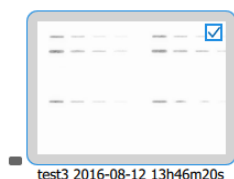

3. In the bottom toolbar, tap View Selected Images to display the selected image in Image View. The image you selected appears full size in Image View.
4. In the Image View bottom toolbar, tap Image Info to open the Image Information box.
5. Under Acquisition Information, obtain the Exposure Time (sec) for the selected image.

Acquisition Information	
Imager	ChemiDoc™ Touch MP
Exposure Time (sec)	120.000 (Signal Accumulation)
Serial Number	KNOWNMP
Software Version	0.9.0.1225
Application	Chemiluminescence
Excitation Source	No Illumination
Emission Filter	647SP Filter
Binning	2x2

6. Close the Image Information box.

To acquire an image using the best SAM exposure time

1. Tap Camera view.
2. Tap Exposure and then tap Manual.
3. Tap Set Manual Exposure Time.

4. Tap the sec (seconds) box to display the on-screen keypad.
5. Enter the Exposure Time value you obtained from the Image Information box.
6. Tap Camera to acquire the image.

Adjusting SAM Acquisition Settings

You can preview SAM images as they are acquired and cancel the acquisition at any time. You can then adjust the SAM settings and start SAM acquisition with the adjusted settings.

When you cancel a SAM acquisition in progress, the imager stops acquiring images and stores the images already acquired in the Gallery. For example, if you set the total number of images to 10 and cancel while image 4 is being acquired, acquisition stops immediately, and images 1–3 are stored in the Gallery.

To cancel SAM acquisition and adjust acquisition settings

1. Carefully preview each image as it is acquired.
2. If an image is not to your liking, tap Cancel.

The software stops the acquisition, stores acquired images in the Gallery, and displays the last image acquired at full size.

3. Go to Camera view and tap Exposure.
4. Change the SAM settings.
5. Tap Camera to start another SAM acquisition.

Acquiring the Image

After you specify the image settings, you can acquire the image. Make sure the transilluminator drawer is pushed in all the way and the door is closed.


Note: If an exclamation point appears on the screen in place of the Camera, a problem is preventing image acquisition. The color of the exclamation point indicates the severity of the problem. An orange exclamation point indicates that an action must be taken before the image can be acquired. A red exclamation point indicates that the camera is disconnected. Log out of Image Lab Touch and restart the imager to correct the error.

To view the error message

- ▶ Tap the exclamation point button in the bottom left corner of the screen.



A screen message explains how to correct the error.

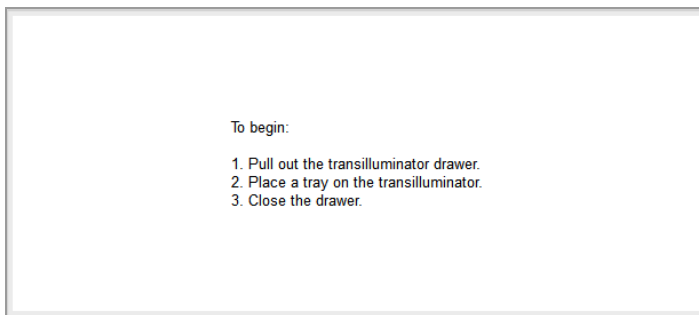

To acquire a gel or blot image

- ▶ Tap Camera.

Acquisition starts. A progress bar monitors the acquisition.

When the image is acquired, it appears in Image View. You can adjust the appearance of the image. You can also print, rename, and delete the image in this view.



(Multichannel) If you cancel the process during multichannel acquisition, the software stops acquiring images and stores the images that have been acquired in the Gallery.

(Stain-Free Gel) Skipping Gel Activation

When you select a gel activation time for a stain-free gel and tap Camera, gel activation starts before acquisition. A preview of the image appears. You can monitor the progress of gel activation and stop it when the gel is properly activated.

To stop gel activation and skip to acquisition

1. Monitor gel activation carefully.
2. When the image is properly activated, tap Skip in the progress box.

The software skips the remaining time for gel activation and acquires the image.

(Stain-Free Gel) Acquiring an Image after Canceling Acquisition

Important: Do not open the door or move the sample.

When you activate a stain-free gel and then cancel image acquisition in progress, you can use the same gel, keep or change exposure settings, and acquire the image again.

To acquire an image after canceling acquisition

1. Tap Exposure.
2. Select No Activation.
3. Set exposure options.
4. Tap Camera to reacquire the image.

(Multichannel) Retaking an Image

You can retake one or more channel images in a multichannel image right after acquisition. You can also delete one or more channels to generate a different multichannel image. The new multichannel image results from the combination of the remaining channel images.

Important: Do not leave the current view, open the door, or log out. Any of these actions makes retaking an image unavailable.

You can change the exposure settings and retake the channel image to achieve different results. You can also change automatic or manual exposure options.


When you retake an image, Image Lab Touch does the following:

- Replaces the channel image you specify with a new image acquired using modified exposure settings. The other channel images are retained.

- Generates a new composite image by combining the channel image you retake with the channel images you keep. (A *composite image* combines the data from the channel images into one image.)
- Stores the original multichannel image and the new multichannel image in the Gallery. (A *multichannel image* consists of a composite image plus the channel images generated during one acquisition.)


To retake a channel image

1. Tap Retake in the main toolbar.

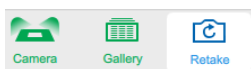

2. Change the exposure settings in one or more channels.


The previous and current settings appear under Exposure in the imaging settings tile.

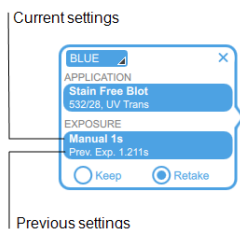

3. (Optional) Select an automatic exposure and change the region of interest.
4. (Optional) To revert back to the previous settings, tap Keep in the imaging settings tile.
5. To retake the image, tap Camera.

To delete a channel and retake a multichannel image

1. Tap Retake in the top toolbar.
2. Tap the X in the upper right corner of the imaging settings tile to delete the channel and the channel image.

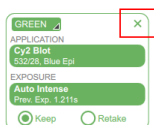

3. To retake the image, tap Camera.

The software generates a new composite and multichannel image.

Merging Images


You can merge images to create multichannel images from up to three single images already acquired. For example, you can merge a colorimetric image of prestained standards and a chemiluminescent image of the same blot into a multichannel image.

Original image data are not affected and can be used for quantitation as well as for determining the molecular weight or size of proteins. You can view merged images in gray scale. File names for merged images use the format specified in User > File Naming Preferences on the main toolbar.

Note: You cannot merge multichannel images. You cannot merge a SAM group with another SAM group or with a single image.


To merge images

1. In the Gallery, select up to three images. In this example, two images are selected.

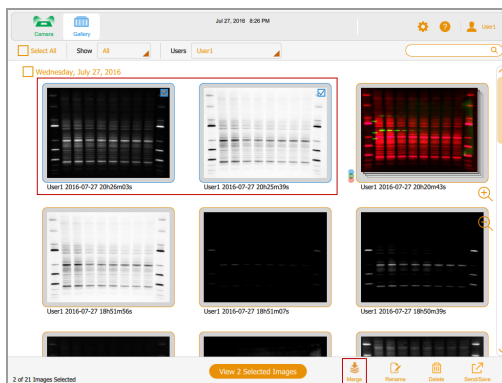


2. On the lower toolbar, tap Merge.

The merged image appears in Image View.

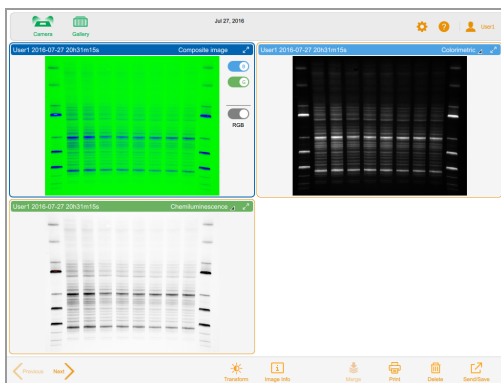

To view merged image information

- On the Gallery lower toolbar, tap Image Info. In the Image Information box, detailed information about each image merged appears in Notes.

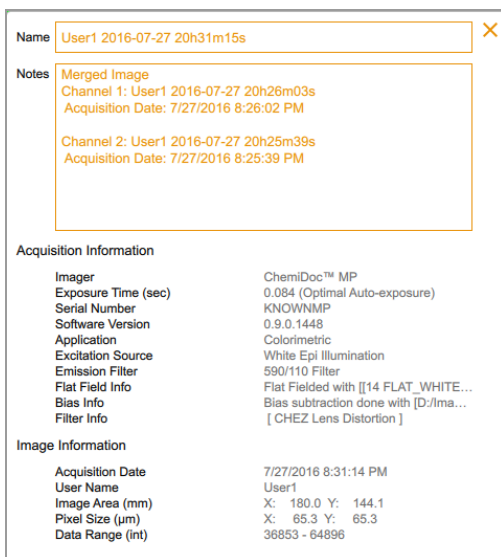

Naming Merged Images

The system names the merged image by combining the names of both images separated by a plus sign.

Example

2015-04-30 17hr 26min 19sec+2015-04-30 17hr 27min 24sec

If the name is already in use in the Gallery, the system generates another name by adding a hyphen and an incremental number (starting at 1) to the end of the combined name. The system continues to increase the added number incrementally (-2, -3, -4 and so on) until the name is unique.

Example


2015-04-30 17hr 26min 19sec+2015-04-30 17hr 27min 24sec-1

The system stores the names of the merged image and the two source images.

To view the names of the source images

- With the merged image open in Image View, tap Image Info in the bottom toolbar.

In the Image Info box, the Notes section displays the names of the images that were merged.

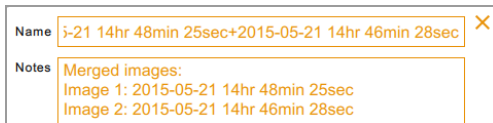

Chapter 7 Excising Bands from Samples

You can excise bands of interest from agarose or acrylamide gels for applications such as mass spectrometry or DNA cloning.

The procedure for excising bands varies depending on the sample tray in use.

Excising Bands on a White or Blue Tray

Working with white and blue trays does not require using the UV shield. However, you must wear yellow XcitaBlue goggles to see bands on a blue tray.



Caution: Sharp cutting tools can easily damage the surface of the trays. To avoid this, place a sheet of clear glass or plastic on the tray before you add the gel sample. Use a *chopping* motion rather than a *sawing* motion.

To excise bands on a white or blue tray

1. Place a gel sample in the center of the tray.
2. In Camera view, tap Turn Transilluminator On.

The transilluminator turns on, illuminating the gel.

Note: The lights turn on only when the sample tray is in place.

3. Excise the bands.
4. When you have finished excising the bands, tap Turn Transilluminator Off.
5. Slide in the transilluminator drawer and close the door.

Excising Bands on a Blot/UV/Stain-Free Tray




WARNING! Transilluminators are powerful sources of UV radiation, which can cause serious damage to unprotected eyes and skin. The accessory UV shield (catalog #1708375) provides some UV protection. However, this shield does not protect others standing in the area around the imager. Before performing band excision, the user and other lab personnel in proximity to the imager must put on protective gear including UV protective safety glasses, a face shield, lab coat, and gloves to ensure that no skin is exposed. A typical and reasonable expectation of use is three operations a day per user for three minutes each. Bystanders without protective gear must stand at least 1.5 meters (five feet) away from the imager and limit their exposure to no longer than one hour per day.


Important: Before excising bands, you must install the UV shield.

To install the UV shield


1. Open the imager front door and pull out the transilluminator drawer.
1. Open the imager front door and pull out the transilluminator drawer.
2. Place a UV tray on the transilluminator drawer.
3. Hold the UV shield by diagonal corners so you can guide it accurately.

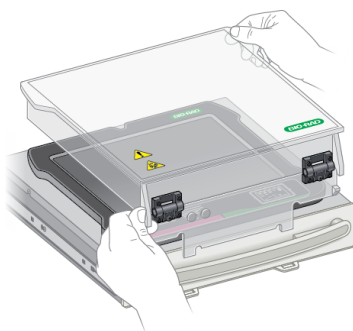


4. Insert the two locking tabs that extend from the front of the UV shield into the two slots on the front of the transilluminator drawer.

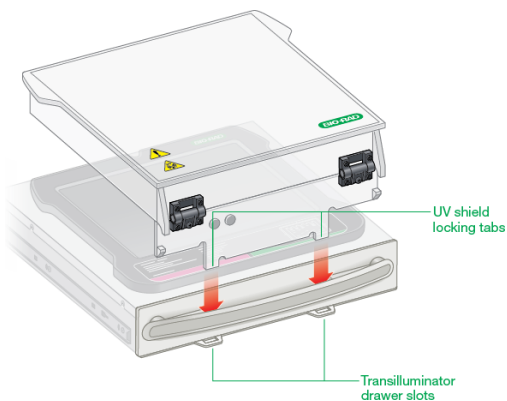

5. Ensure that the two wedges on the front of the UV shield snap into position on the transilluminator handle.

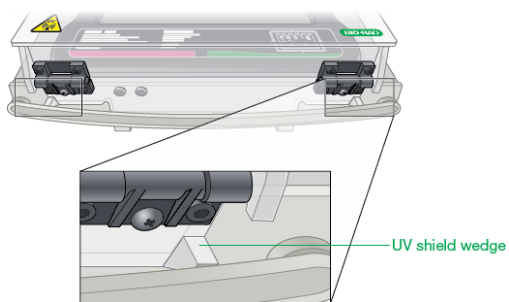

The installed UV shield should look like this:

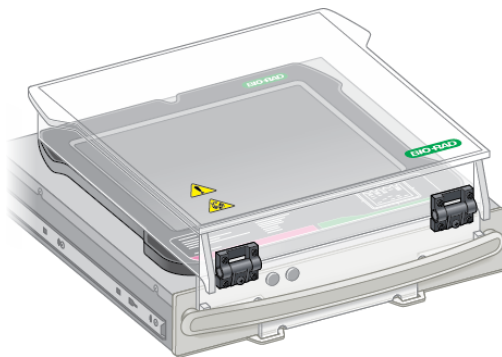

Important: Before you begin excising bands, you must put on the required protective gear and ensure that the UV shield is installed.

The UV lights turn off after 15 minutes of continuous use. To turn the UV lights back on, tap Turn Transilluminator On.

To excise bands on a Blot/UV/Stain-Free tray

1. To avoid damaging the surface of the sample tray, place a sheet of clear glass or plastic on the tray before you add the gel sample.
2. Place a gel sample in the center of the tray.
3. In Camera view, tap Turn Transilluminator On.

Turn Transilluminator On changes from green to blue and the UV lights turn on, illuminating the gel.


Note: The UV lights turn on only when both the sample tray and the UV shield are in place. If the lights do not turn on, verify that the UV shield is installed correctly.

4. Raise the UV shield no more than is necessary to work with the sample.



Caution: Keep the UV shield open for as little time as possible.

5. Reach around the sides of the shield to excise the bands.

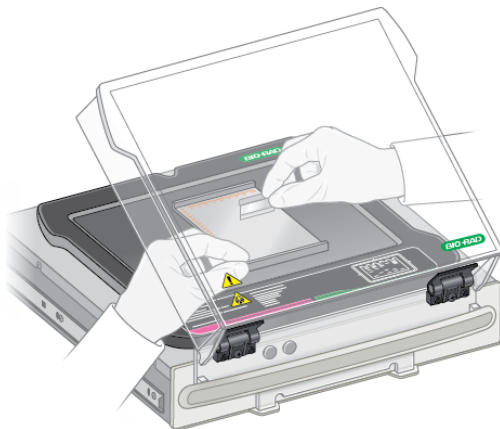

Caution: Sharp cutting tools can easily damage the surface of the tray. Use a *chopping* motion rather than a *sawing* motion.

6. When you finish excising bands, tap Turn Transilluminator Off to turn off the UV lamps.
7. Close the UV shield.
8. Remove the UV shield, remove the sample tray, slide in the transilluminator drawer, and close the door.

Chapter 8 Viewing Images

You can view images in Camera view or in the Gallery. The Gallery displays thumbnails of all images you acquire, in order by the date of acquisition, most recent acquisition date shown first.


Note: Administrators can view all users' images at once or view an individual user's images.

Accessing the Gallery

To access the Gallery

- Tap Gallery in the main toolbar.

A Gallery similar to the following appears.

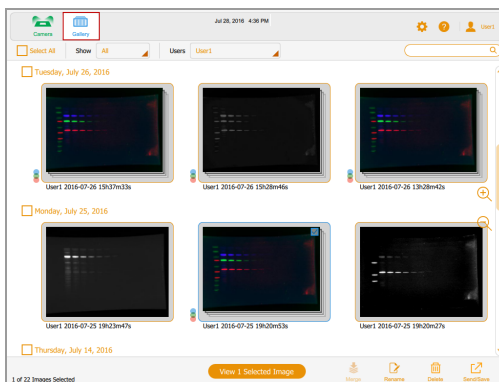

You can change the size of the thumbnails in the Gallery by tapping the plus and minus zoom icons on the right side of the Gallery screen.

The image name appears below each thumbnail image. By default, each name consists of the user name and the date and time of acquisition, for example: User1 2016-07-16 15hr7min33sec.scn.

You can change the file name in the Image Information box. You can also change the image name format. See [Changing the File Naming Format on page 132](#).

To open an image in the Gallery

- Double-tap a thumbnail image.

The image opens full size in Image View.



Selecting and Opening Images in the Gallery

To select an image

- ▶ Tap a thumbnail.


The selected image thumbnail appears outlined in blue and a check box appears in the upper right corner.

To open the selected image

- ▶ Tap the selected image to open it full size in Image View.

To open 1–4 images

- ▶ Select 1–4 images and tap View x Selected Images in the bottom toolbar (where x is the number of images selected).

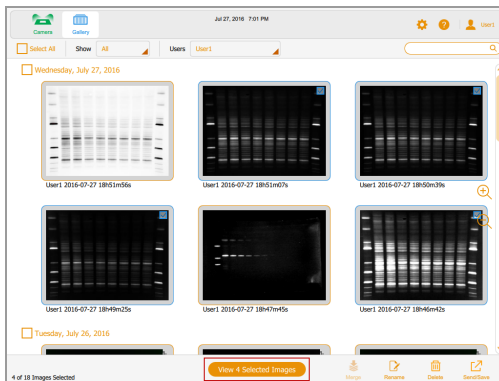


The selected images appear outlined in blue.

To select all images you acquired on a single date


- ▶ Tap a date checkbox.

All images you acquired on the selected date appear outlined in blue.

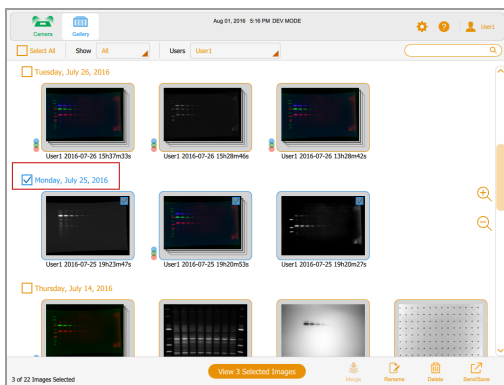

To select all of your images


► Tap Select All.

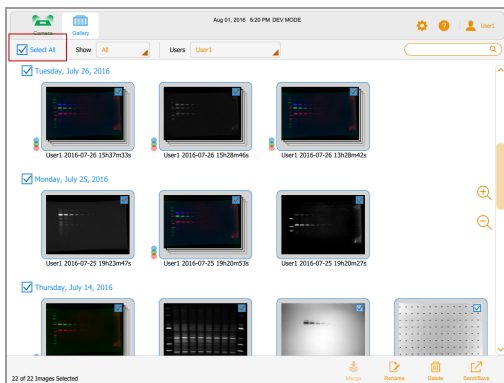

All images are selected and appear outlined in blue.

Filtering Gallery Images by Date

By default, the Gallery shows all images you acquired starting with the last image you acquired on the date you specify. When you acquire another image, the Gallery displays that image and all other images in the Gallery.

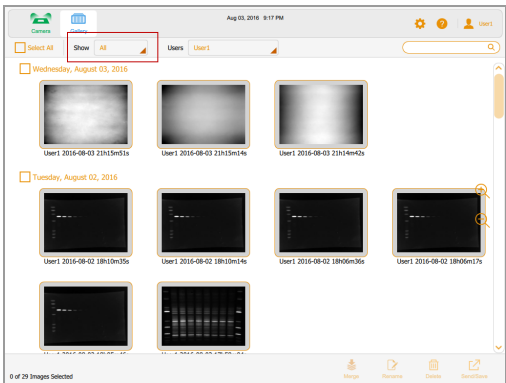


You can filter the display by date to focus on images of immediate interest.

To filter images by date

1. Tap Show All in the image selection toolbar.

A dropdown list of date choices appears.

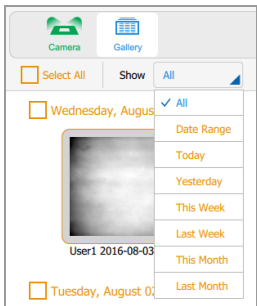

2. Tap one of the following.

- **Date Range** — displays images acquired within a range of dates you specify. Tap the calendar boxes to specify dates in any order.



- **Today** — displays images you acquired today starting at 12:00:00 AM and ending at 11:59:59 PM
- **Yesterday** — displays images you acquired yesterday starting at 12:00:00 AM and ending at 11:59:59 PM.
- **This Week** — displays images you acquired this week starting at, for example, 12:00:00 AM Sunday and ending at 11:59:59 PM the following Saturday.
- **Last Week** — displays images you acquired starting at 12:00:00 AM Sunday last week and ending at 11:59:59 PM the following Saturday.
- **This Month** — displays images you acquired this month starting at 12:00:00 AM of the first calendar day of the month and ending at 11:59:59 PM of the last calendar day of the month. For example, the month of January starts at 12:00:00 AM January 1 and ends at 11:59:59 PM January 31.
- **Last Month** — displays images you acquired starting at 12:00:00 AM of the first calendar day of last month and ending at 11:59:59 PM of the last calendar day of last month.

Checking Intensity Values

In Image View, you can display the average intensity value and the percentage of saturation for any spot in an image. A 3 x 3-pixel region of the image is averaged, except at the edge or corner, where a 3 x 2 or 2 x 2-pixel region is averaged.

When viewing a multichannel image, you can check intensity values in any of the channel views or in the composite view.


Note: Hiding a channel in a composite image also hides the intensity values for that channel.

You can use the saturation percentage to estimate how much to increase exposure time to maximize the dynamic range of the image. For example, if the spot is 50% saturated, doubling the exposure time results in about 100% saturation. Viewing intensity values is also useful when comparing images.


To view intensity values in an image

1. In Image View, display an image.
2. Tap any spot in the image.

Intensity values for the selected spot appear.

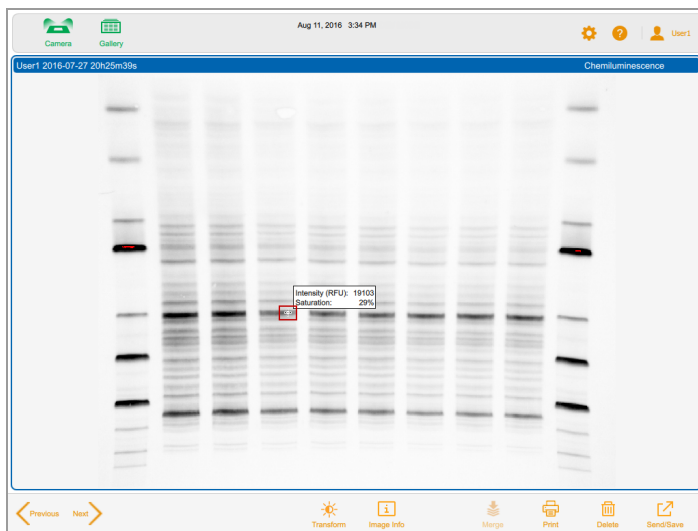


Crosshairs mark the pixel area being evaluated. Intensity values fade out after a few moments.

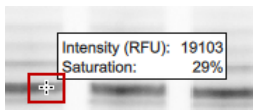

To view intensity values in a composite image

1. In Image View, display a multichannel image.
2. Tap any spot in the composite image.

Color coded intensity values appear for each channel in the composite image.

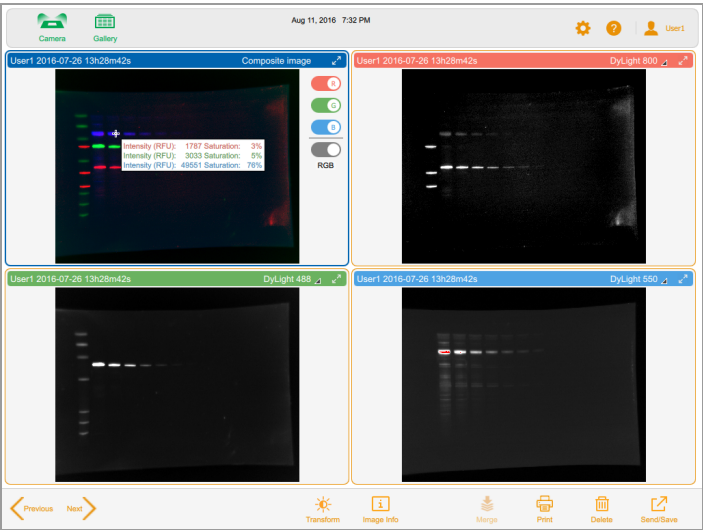


Crosshairs mark the pixel area being evaluated. Intensity values fade out after a few moments.



Viewing Multichannel Images in Image View

After a multichannel image is acquired, the composite image and its individual channel images appear in Image View. A blue border surrounds the selected image.

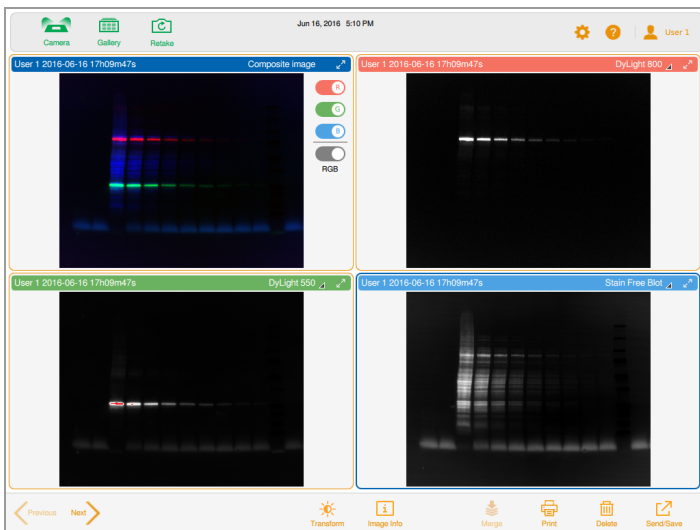


Images you open in the Gallery also appear in Image View. You can select and open up to four images at a time in the Gallery, all of which appear together in Image View.

To view more than one multichannel image at a time


1. In the Gallery, select multichannel images you want to view.
2. In the lower toolbar, tap View x Selected Images, where x is the number of images you selected.

The multichannel images you selected appear in Image View.

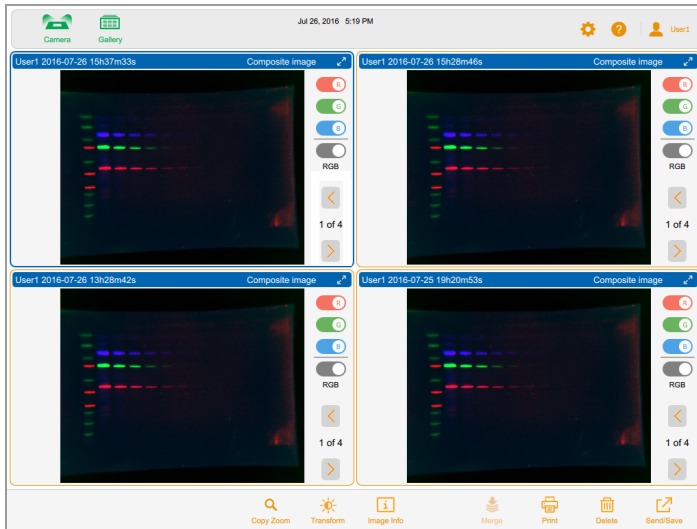


To view a single image in a multichannel image

- Tap the Previous and Next arrows below the RGB settings.

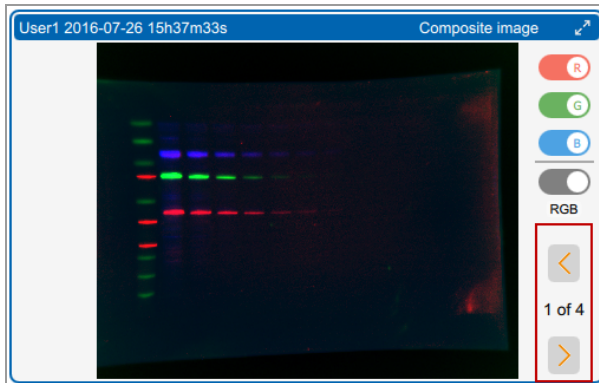

To view an image full size


- ▶ Tap the expand arrows in the upper right corner of the image.

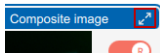

Viewing a Composite or Merged Image in Gray Scale

In Image View, you can view the composite or merged image in gray scale. Channel colors of individual images are unaffected by the gray scale setting. You can also export or print a composite or merged image in gray scale. In the Gallery, the thumbnail of the image also appears in gray scale.


By default, a composite or merged image appears in Image View in RGB, the combined color settings of its channels.

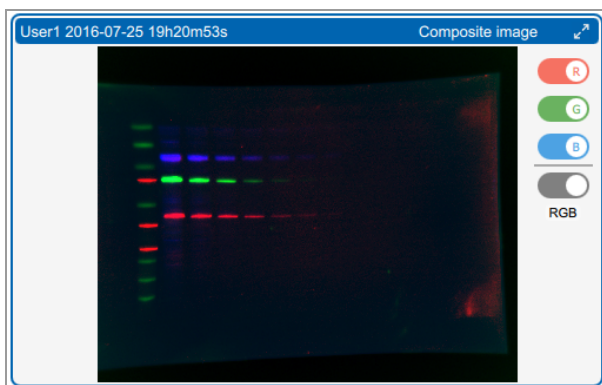

To view the image in gray scale

- ▶ Tap the RGB toggle below the channel colors.

The toggle name changes to Gray and the image appears in gray scale.

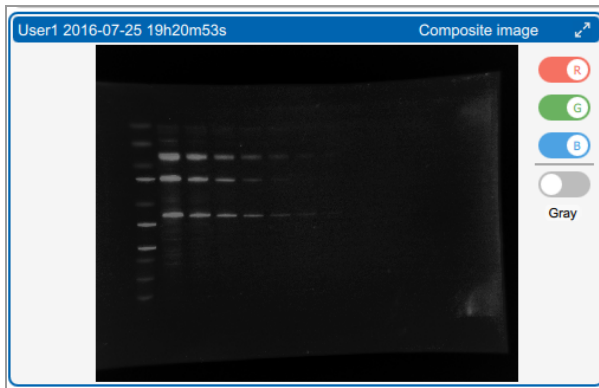

Comparing Images


In the Gallery you can open and compare up to four images of a gel or blot at a time.

To open multiple images

1. Select 1–4 images in the Gallery.
2. In the bottom toolbar, tap View x Selected Images (where x is the number of images selected).

The images you select in the Gallery open in Image View. By default, the top left image is selected.

Chapter 8 Viewing Images

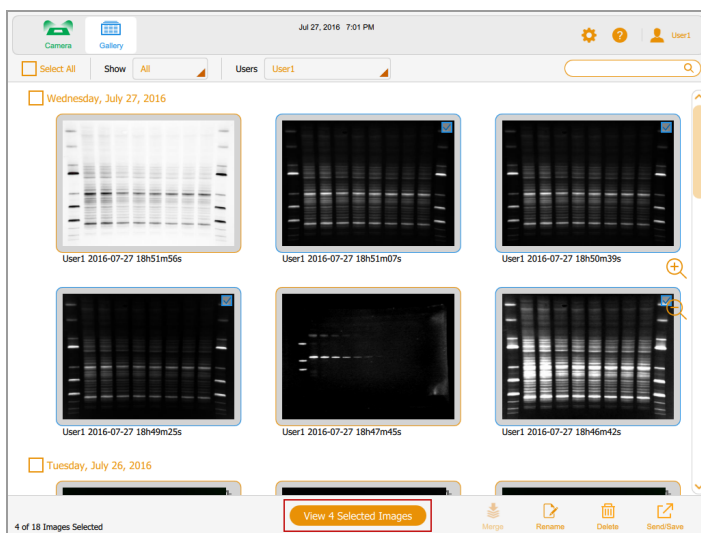


Copying Image Settings

It is easier to compare images that have the same transform and zoom settings. You can copy these settings from one image to the others you selected.

To copy zoom settings from one image to the others

1. Select the image with the zoom settings you want to copy.

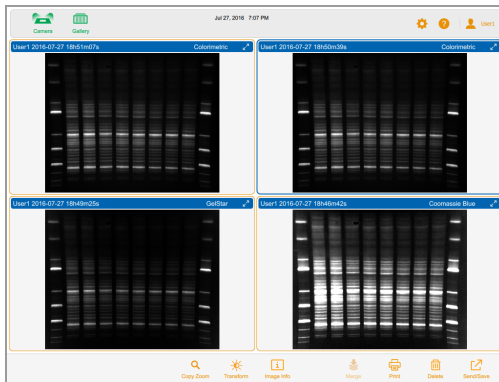

2. In the bottom toolbar, tap Copy Zoom.




The pan and zoom settings of the selected image are copied to the other images.

To copy transform settings from one image to the others

1. Select the image with the transform settings you want to copy.
2. In the bottom toolbar, tap Transform.



The transform settings for the selected image appear.

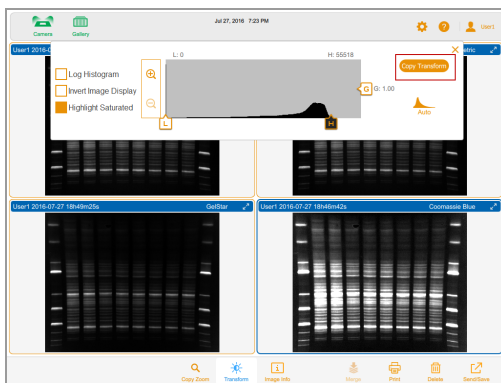

3. Tap Copy Transform in the transform settings box to copy the transform settings to the other images.

Viewing and Editing Image Information

The Image Information box provides information such as the image name, the acquisition's exposure duration and date, and the application type. You can change the image name in the Name box. You can also add information about the image in the Notes box.

To view information about the image

- In Image View, tap Image Info in the bottom toolbar.

The Image Information box appears.

Name

User1 2016-08-09 19h32m45s

×

Notes

Acquisition Information

Imager	ChemIDoc™ MP
Exposure Time (sec)	0.383 (Optimal Auto-exposure)
Flat Field	Applied (Blue Epi)
Serial Number	KNOWNMP
Software Version	0.9.0.1448
Application	Fluorescein
Excitation Source	Blue Epi Illumination
Emission Filter	532/28 Filter

Image Information

Acquisition Date	8/9/2016 7:32:45 PM
User Name	User1
Image Area (mm)	X: 180.0 Y: 144.1
Pixel Size (µm)	X: 65.3 Y: 65.3
Data Range (int)	40895 - 65535

To edit image information

1. Tap inside the Name or Notes box.
2. Make the text changes you want.
3. Tap OK on the on-screen keyboard to save your changes.


Viewing SAM Images

When SAM images are acquired, the last image acquired appears full size in Image View. You can view all the images in a SAM acquisition in the Gallery, where you can also rename, delete, or export the SAM group.


To select a SAM acquisition

1. Tap Gallery.


The Gallery displays the stacked SAM images in a thumbnail. The file name appears below the thumbnail. A stack icon in the bottom left corner identifies the thumbnail as a SAM acquisition.

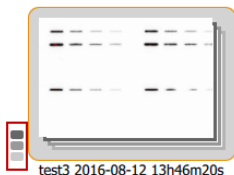

2. Tap the SAM stack thumbnail.

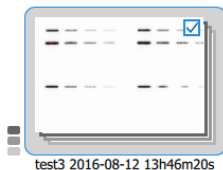


The bottom toolbar displays the number of SAM images in the acquisition:

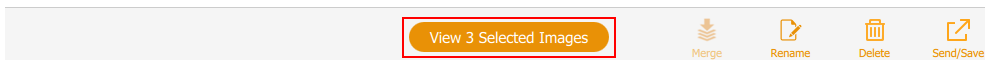


To view all images in a SAM acquisition

1. In the Gallery, tap the stack icon.

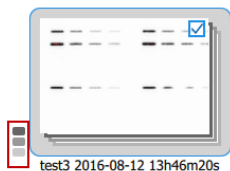

A thumbnail of each Image in the SAM acquisition appears.

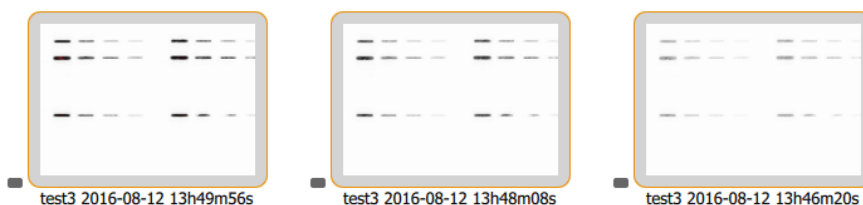

To view a SAM image in Image View

- Double-tap a SAM image thumbnail.

The image you selected appears full size in Image View.


To view up to four SAM images in Image View

- Tap up to four SAM image thumbnails and then tap View x Selected Images in the bottom toolbar, where x equals the number of images you selected.

The selected images appear in Image View.

To restack SAM images

- ▶ Tap the SAM icon on any thumbnail in the SAM acquisition.

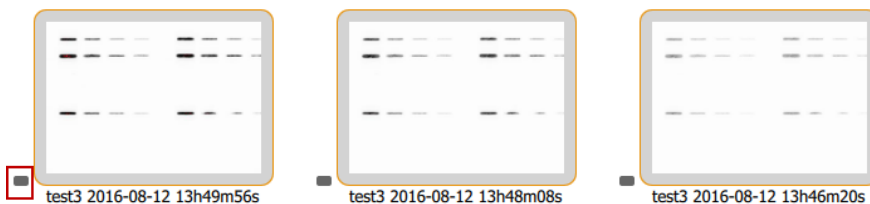

To clear a SAM image selection

- ▶ Tap inside a SAM image thumbnail.

To clear all image selections and restack the SAM images

- ▶ Clear the checkbox in any thumbnail and tap its stack icon.

Chapter 9 Managing Images

Adding Notes to an Image File

To add information to an image file

1. In Image View, tap Image Info in the bottom toolbar.

The Image Information box appears.

The screenshot shows a dialog box titled 'Image Information'. It has a 'Name' field at the top with the text 'User1 2016-07-08 14h56m58s' and a close button (X) to its right. Below the name field is a large 'Notes' text area. Underneath the notes area is a section titled 'Acquisition Information' which contains a list of technical details in two columns. At the bottom is a section titled 'Image Information' which contains a list of acquisition parameters in two columns.

Acquisition Information	
Imager	ChemIDoc™ MP
Exposure Time (sec)	0.042 (Optimal Auto-exposure)
Serial Number	KNOWNMP
Software Version	0.9.0.1418
Application	Colorimetric
Excitation Source	White Epi Illumination
Emission Filter	590/110 Filter
Flat Field Info	Flat Fielded with [14 FLAT_WHITE...
Bias Info	Bias subtraction done with [Dz/Ima...
Filter Info	[CHEZ Lens Distortion]


Image Information	
Acquisition Date	7/8/2016 2:56:56 PM
User Name	User1
Image Area (mm)	X: 180.0 Y: 144.1
Pixel Size (µm)	X: 65.3 Y: 65.3
Data Range (int)	32205 - 64603

2. Tap in the Notes box.
3. Enter your notes using the on-screen keyboard.
4. Tap OK to save the notes and close the keyboard.

For more information, see [Viewing and Editing Image Information on page 126](#).

Changing the File Naming Format

By default, the file name consists of the elements user name, date, and time. In the File Naming Preference dialog box, you can change the file naming format at any time.

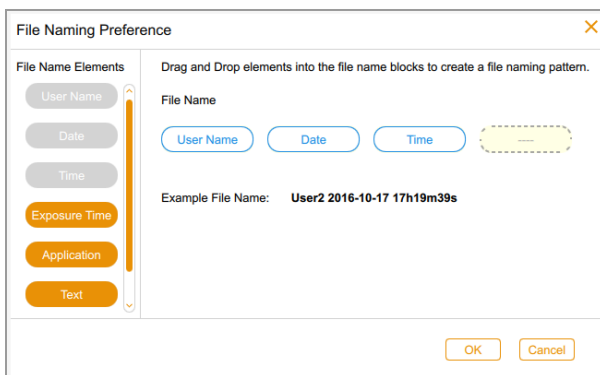


- You can remove an element and replace it with another element.
- You can include the Text element, which holds up to 32 alphanumeric characters.
- You can reorder the elements.

To change the file naming format, you select up to four of the seven elements in the File Name Elements List and arrange them in the order you prefer.

Tip: The File Name Elements list includes a combination element, App/Exposure, which is useful when you want to include more than four kinds of information in the format.

To access the File Naming Preference dialog box

- Tap User Settings, and then tap File Naming.

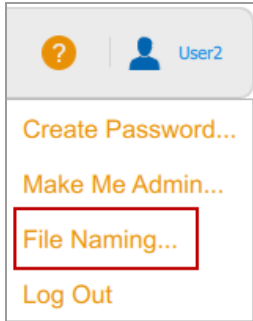

The File Naming Preference dialog box appears.

To change the file naming format

1. In the the File Naming Preference Dialog box, drag an element you want to exclude from the format onto the File Name Elements list.
2. Scroll down the File Name Elements list, choose an element you want to include, and drag it into position in the file name.

An example of the file name in progress appears as you select each element.

3. Repeat steps 1 and 2 until the file name format includes all the elements you want.
4. (Optional) Rearrange the elements to your liking.
5. Tap OK to save the format.

The software names all image you acquire using the new file naming format until you change it.

Renaming Images

An image name can consist of up to 190 alphanumeric characters. Spaces are permitted except at the beginning and end.

Note: The following characters cannot be used: \ / : * ? " < > |

To rename an image in the Gallery

1. Select an image.
2. Tap Rename in the bottom toolbar.
3. In the Rename dialog box, enter a name for the image.
4. Tap OK.

To rename multiple images in the Gallery

1. Select more than one image.
2. Tap Rename.
3. In the Rename dialog box, enter a name for the first image.
4. Tap OK.

The system labels the remaining images with the same name as the first image plus a consecutive number. For example: ChemiBlot_1, ChemiBlot_2, and so on.

To rename an image in Image View

1. In Image View, tap Image Info.
2. In the Image Information box, tap in the Name box to select the image name.
3. Type a new name.
4. Tap OK to save the image name.

Renaming SAM Images

You can rename a SAM image or a SAM group. You can rename a combination of single SAM images, SAM groups, and other images. Image and group names must be unique.

When you rename a SAM group each image in the group is also renamed. An underscore and a sequential number are added to each image name to render it unique. The oldest SAM group or image is assigned the first sequential number. The remaining images are assigned sequential numbers, oldest image first.

Example

You select a SAM group, SAM 8497, and another image in the Gallery.

You rename SAM 8497 to SAM 9001.

The software adds the suffix _1 to the renamed SAM group. The thumbnail for the SAM group now reads SAM 9001_1.


Each image in group SAM 9001_1 is also renamed. The first image acquired is named SAM 9001_2. The remaining images (including the image that is not in the SAM group) are renamed SAM 9001_3, SAM 9001_4, SAM 9001_5, and so on, in the order in which they were acquired. If the image outside the SAM group was acquired before the acquisition of the SAM group, that image is renamed SAM 9001_1. In this case, the SAM group is renamed SAM 9001_2. The remaining images are renamed SAM 9001_3, and so on, in the order in which they were acquired.

To rename a SAM image

1. Tap the SAM stack thumbnail.
2. Select a SAM image.
3. Tap Rename in the bottom toolbar.
4. Enter a name in the Rename Image dialog box and then tap OK.

To rename all images in a SAM group

1. Restack the SAM images.
2. Tap the SAM stack icon.
3. Tap Rename.
4. Enter a name in the Rename Image dialog box and tap OK.

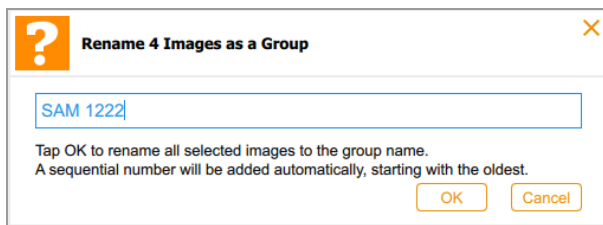

Printing an Image

You can print one image at a time on the Mitsubishi P95 printer only. The printer produces a black and white image in landscape orientation. The printout includes the user name, acquisition date and time, and the application used.

To print an image

1. Ensure that the printer is connected to the imager and is turned on.
2. Do one of the following:
 - In Image view, display the image you want to print.
 - In the Gallery, select the image you want to print.
3. Tap Print.


For more information about the printer, see the operation manual for your printer.

Searching for Images in the Gallery

You can search for images of interest in the Gallery. Search text is not case sensitive. You can enter any part of the date or name. You can also search by application name or type.

To search for an image

1. Tap the search box to display the keyboard.

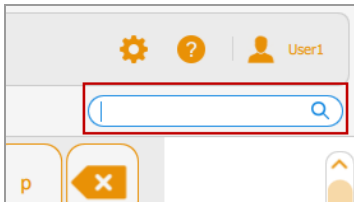

2. Enter search text.

As you enter search text, images that match the search text appear in the Gallery.

3. Tap Done.

To return to a thumbnail view of all your images

- Clear the search box.

Deleting Images

In Image View, you can delete any open image. In the Gallery, you can delete multiple images at the same time.

To delete images in the Gallery

1. Do one of the following to select images to delete:
 - Tap one or more thumbnails.
 - Tap the checkbox for one or more days to select all thumbnails for those days.

- Tap Select All to select all thumbnails in the Gallery.
2. Tap Delete and then tap OK.

To delete an image in Image View

1. Do one of the following:
 - If one image is open in Image View, tap Delete.
 - If more than one image appears in Image View, select the image you want to delete.
2. Tap Delete and then tap OK.

Deleting SAM Images

You can delete SAM images in the Gallery or in Image View.

In the Gallery, you can delete

- A SAM image or a SAM acquisition
- A combination of single SAM images, SAM acquisitions, and other images

In Image View, you can delete

- A full size image
- An image you select from a view of up to four images

To delete images from a SAM acquisition in the Gallery

1. Tap the stack icon next to a SAM stack thumbnail.
2. Tap the thumbnail of one or more SAM images you want to delete.
3. Tap Delete in the bottom toolbar.
4. To confirm deleting the images permanently, tap OK.

To delete a SAM acquisition in the Gallery

1. Restack the SAM images.
2. Tap inside the SAM stack thumbnail.
3. Tap Delete in the bottom toolbar.
4. To confirm deleting the images permanently, tap OK.

To delete a combination of SAM images, other images, and SAM acquisitions in the Gallery

1. Tap the stack icon next to each SAM acquisition that contains images you want to delete.
2. Tap the SAM image thumbnails, non-SAM images, and SAM acquisitions you want to delete.
3. Tap Delete.
4. To confirm deleting the images permanently, tap OK.


To delete a SAM image in Image View

1. Do one of the following:
 - If one SAM image appears in Image View, tap Delete.
 - If more than one image appears in Image View, select the image you want to delete and tap Delete.
2. To confirm deleting the image permanently, tap OK.

Chapter 10 Exporting Images

After acquiring images using Image Lab Touch software, you can import your images into Image Lab. To do so, first export the acquired images to a USB flash drive or a shared folder. Then copy the exported images to a computer running Image Lab software.

You can export a SAM group or individual SAM images. You can also export a combination of single SAM images, SAM groups, and other images. SAM images and groups are stored in a folder inside the top-level time-stamped folder:

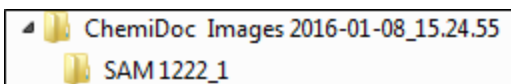

You can export images from the imaging system to one of the following:

- Flash drive or external hard drive
- Shared folder on a network drive or stand-alone computer

Image Lab Touch exports the images to a folder at the top level of the destination location. The folder is named *ChemiDoc Images* plus a time stamp, for example, ChemiDoc Images 2017-04-29_19.08.51. A folder with a time stamp is created each time you export images.

For More Information

For more information about image analysis, see the Image Lab Software User Guide.

Preparing to Export Images

Before you can export images to a stand-alone computer, a shared folder must be present on the computer, and you must provide the UNC path to that folder.

The UNC (universal naming convention) is a standard for identifying servers and computers on a shared network. It is the full network pathname of a folder, including the network server name. This path is the same as a Windows folder path except that it uses double backslashes (\\) to precede the name of the computer.

Note: The procedures for determining the UNC path and creating a shared folder on it are based on recent versions of Windows and Mac operating systems. If your computer is running on a different operating system version, you might notice differences in the user interface.

Creating a Shared Folder on a Windows Computer


This section explains how to create a shared folder on a Windows computer and how to determine its UNC.

Note: In order to export images to a Windows computer, ensure that the imaging system is connected to the stand-alone computer or to the network port by Ethernet cable.

To create a shared folder on a Windows computer

1. Log in as the Windows Administrator for the computer.
2. Create a destination folder.
3. Right-click the folder and select Properties in the menu that appears.
4. Click the Sharing tab in the Properties dialog box.


5. Under Network File and Folder Sharing, click Share.

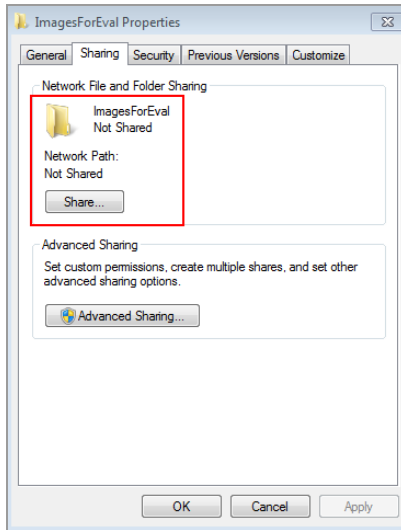

6. Choose users with whom you want to share the destination folder.
7. Click Share and then click Done.
8. Set Read/Write permissions for each user.
9. Click Close.

To determine the UNC path to the shared folder

- On the Sharing tab, in the Network Files and Folder Sharing section the UNC path appears under Network Path.

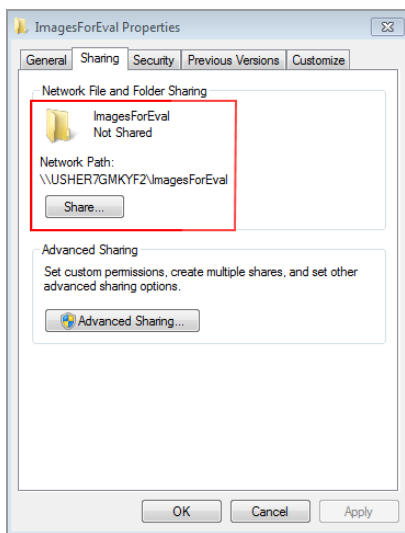

Write down this path. You will enter this address when you export an image.

Creating a Shared Folder on a Mac Computer

This section explains how to create a shared folder on a Mac computer and how to determine its UNC.

Tip: The UNC for a Mac computer includes its Apple Filing Protocol (AFP) or Server Message Block (SMB).


Note: In order to export images to a Mac computer, ensure that the imaging system is connected to the stand-alone computer or to the network port by Ethernet cable.

To create a shared folder on a Mac computer

1. Log in as the Administrator.
2. Create a destination folder.
3. Click the folder and choose Get Info in the menu that appears.


The *<folder name>* Info dialog box appears.

4. Select Shared Folder.

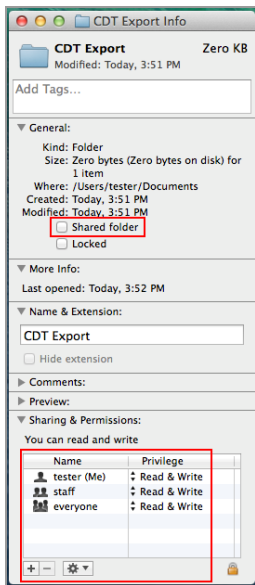

5. In the Sharing & Permissions box, verify that the users with whom you want to share the folder have permission to read and write to this folder.

6. On the Apple menu, open System Preferences and select File Sharing in the Sharing dialog box.

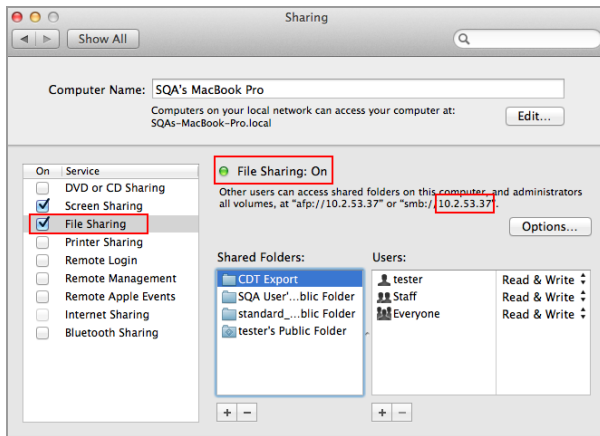

7. In the Service pane, select File Sharing.
8. Under File Sharing: On, locate the SMB address and write it down.
You will enter the SMB address when you export an image.
9. Click Options.
10. In the dialog box that appears, select Share files and folders using SMB.

You can select both SMB and AFP. By default, OS X Mavericks and later systems automatically enable SMB and AFP for compatibility with Windows computers, Macs using Mavericks and Yosemite, and Macs using older versions of OS X.



11. Click Done.

Determining the UNC of a Previously Shared Folder

In order to export images to a computer, you must provide the UNC path to the folder. This section explains how to determine the UNC of a folder that has already been shared on Windows or Mac computer.

Determining the UNC Path to a Network Folder on a Windows Computer

If the folder is located on a mapped network drive, you will need to determine the path to that drive as well. Mapped network drives appear in Windows Explorer as simulated drives. This is indicated by a drive letter, for example (H:). This section also explains how to determine the path to a mapped network drive on a Windows computer.

To determine the UNC path to network folder on a Windows computer

1. Log in as the Windows Administrator for the computer.
2. In the Navigation pane, click the Network icon and navigate to and select the destination folder.

The address bar displays the network path to the destination folder, beginning with the word Network, for example:

Network > USHER7GMKYF2 > ImagesForEval

3. Click the address bar. Then UNC path appears in the address bar, for example:

\\ ImagesForEval

4. Write down this path. You will enter this address when you export an image.

To determine the path to a mapped drive

1. Log in as the Windows Administrator for the computer.
2. In the Navigation pane, click the network folder that contains the destination folder.

The address bar displays the mapped network path to the destination folder, beginning with the word Computer, for example:

Computer > jnorby (\\usherfs\users) (H:) > ImagesForEval

3. To construct the UNC path, follow this rule:

\\<shared_folder_name\mapped_foldername\distination folder

In this example

- \\usherfs\users — is the shared folder location
- jnorby — is the mapped folder name
- ImagesForEval — is the distination folder

For example:

\\usherfs\users\jnorby\ImagesForEval


Determining the UNC Path to a Network Folder on a Mac Computer

Tip: The UNC for a Mac computer includes its Apple Filing Protocol (AFP) or Server Message Block (SMB).

Note: In order to export images to a Mac computer, ensure that the imaging system is connected to the computer or to the network port by Ethernet cable.

To determine the UNC path to a network folder on a Mac computer

1. Log in as the Administrator.
2. On the Apple menu, open System Preferences and select File Sharing in the Sharing dialog box.

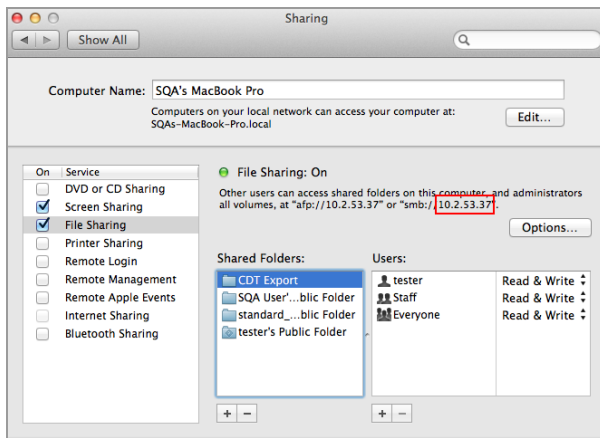

3. Note the AFP or SMB path that appears below File Sharing: On. In this image, the SMB is 10.2.53.37.

This is the computer's IP address, which is used to determine the computer name in the UNC path.

4. To construct the UNC path, follow this rule:

`\\SMB #\Users\<user_name>\<folder_name>`

In this example

- 10.2.53.57 — is the SMB
- Tester — might be the user name
- CDT Export — is the folder name


For example:

`\\10.2.53.57\Users\Tester\CDT Export`

Export Formats

Image Lab Touch exports the image at full size. Zoom and panning views are ignored. The software appends the application name in parentheses to the end of the name of each file saved in .jpg, .tif, or .raw16.tif format. This information is especially helpful for identifying multichannel images.

Example

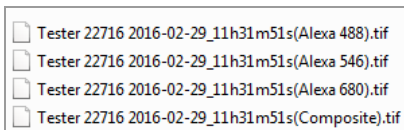

You can export images in the file formats listed on the next page.

Export Option	File Type	Description
Image Lab	.scn or .mscn	<p>Image Lab Touch exports a single-channel image in .scn format or a multichannel image in .mscn format.</p> <p>To work with images in Image Lab software, use these formats.</p> <p>Tip: Because important meta data is exported with your image when exported in these formats,Bio-Rad strongly recommends that you use the .scn or .mscn format.</p>
Publication	.tif or .jpg	<p>Image Lab Touch exports an image in .tif or .jpg format at 300 DPI, which is suitable for publication. You can also insert an image in .tif or .jpg format into presentation software such as Keynote or PowerPoint.</p> <p>Tip: Additional publication options are available in Image Lab software. For example, in Image Lab you can export an image at up to 1200 DPI. To set these publication options for the image, export the image in the .scn or .mscn format, open the file in Image Lab, set the options, and export the image in .tif or .jpg.</p> <p>Note: If you remove a channel from a composite image in a multichannel acquisition and select the .tif or .jpg export option, Image Lab Touch exports the composite image with the remaining channels only.</p>

Export Option	File Type	Description
Analysis	.raw16.tif	<p>Image Lab Touch exports 16-bit .tif format, which retains original pixel intensity data from the image acquisition. This raw image data enhances quantitative analysis. You can also use this format to import the image into image editing software that supports the 16-bit .tif format.</p> <p>Note: You cannot export a composite image from a multichannel acquisition in .raw16.tif format.</p>

Exporting Images


Before you can export images to a network or to a stand-alone computer, a shared folder must be present on the destination computer or network, and you must know the UNC path to the shared folder.

You can export images from the Gallery or from Image View.

To export images


1. Do one of the following:

- Select one or more images in the Gallery to export.

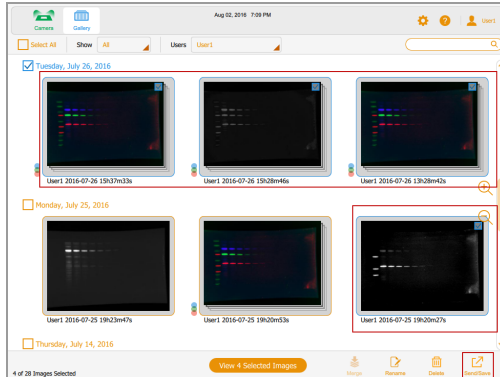

- Display up to four images in Image View and select an image to export.


Chapter 10 Exporting Images

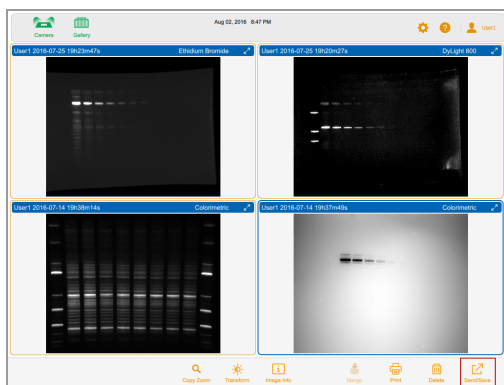

2. Tap Send/Save on the bottom toolbar.


Network and export format settings appear in the lower right corner of the screen.

3. Under Export Options, choose the file formats in which you want to export the images.

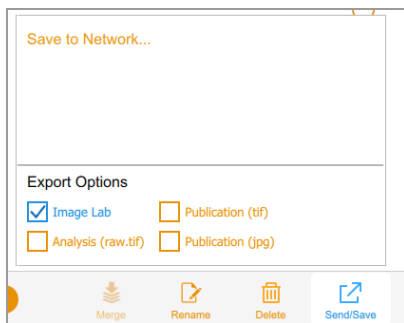

4. In the Network list, select a network.

A progress box monitors the export.

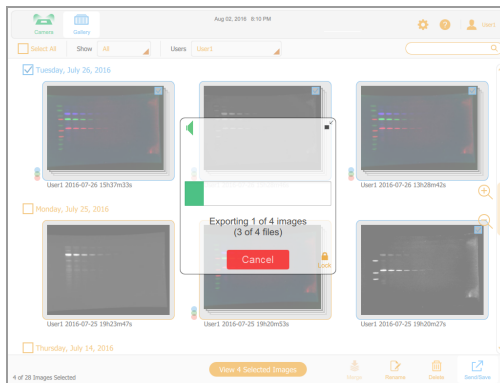

5. When the export finishes, tap OK.

Exporting to a USB Flash Drive or External Hard Drive

The flash drive or external hard drive must meet the following requirements:

- Supports USB 2.0 or above
- Is formatted with FAT or NTFS format
- Has no encryption software or other software add-ons on the drive

Note: USB drives formatted as FAT export files much more quickly than USB drives formatted as NTFS. If your USB export takes too long, reformat the USB drive as FAT.

To export images in the Gallery to a USB flash drive or external hard drive

1. Insert a USB flash drive in or connect an external hard drive to the imager USB port.
2. Do one of the following:
 - Select the images you want to export in the Gallery.
 - Go to Image View by acquiring an image or opening an image in the Gallery.
3. If more than one image is open in Image View, select the images to export.
4. Tap Send/Save on the bottom toolbar.
5. Select export options.
6. Tap Save to USB Drive.

A progress box monitors the image export.

7. When the export finishes, tap OK.

To export images in Image View to a USB flash drive or external hard drive

1. Insert a USB flash drive in or connect an external hard drive to the imaging system USB port.

2. Do one of the following:
 - If the image is displayed in full view, tap Send/Save on the bottom toolbar.
 - If more than one image is displayed, select an image and tap Send/Save on the bottom toolbar.
3. Select export options.
4. Tap Save to USB Drive.

A progress box monitors the image export.
5. When the export finishes, tap OK.

Exporting to a Shared Folder for the First Time

Note: The following characters are not supported in paths to shared folders: * ? " < > | ;

When you enter a path that includes an unsupported character, an error message appears. Folder names that include unsupported characters cannot be selected in browse results. You can rename the shared folder to exclude the unsupported character or choose another folder that does not include one of these characters in its name.


Before exporting images to a shared folder on a network drive or a stand-alone computer for the first time, obtain the following information:

- The UNC path to the shared folder on the network drive or stand-alone computer
- Log-in credentials to the shared folder

To export images to a shared folder

1. Verify that the imager is connected via an Ethernet cable to a network outlet or to a stand-alone computer.

2. Do one of the following:
 - Select the images you want to export in the Gallery.
 - Go to Image View by acquiring an image or by opening an image in the Gallery.
3. In the bottom toolbar, tap Send/Save.
4. Select export options.

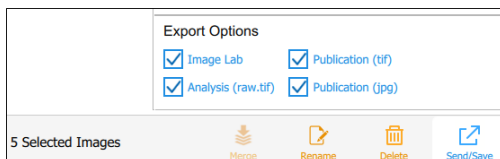

5. Tap Save to Network.
The Save to Network dialog box appears.
6. Type the pathname to the shared folder on the network drive or stand-alone computer.

Enter the network drive server name or IP address.

Server name pathname format: *\\ServerName\Share1\Path\To\Folder*

Example: **\\BioServer1\Share1\Images\Experiment1**

IP address pathname format: *\\IPAddress\Share1\Path\To\Folder*

Example: **\\10.1.52.162\Share1\Images\Experiment1**

7. Tap Save.
8. If a log-in dialog box appears, enter the log-in credentials to the shared folder. For a network drive, enter the network domain name, your user name on the domain, and your password in this form:

domain_name\username\password

Example: **Global.xyz.com\jdoe**

9. Tap OK.

A progress bar monitors the image export. When the export finishes, tap OK.

Exporting to a Shared Folder on a Network or Computer

Note: The following characters are not supported in paths to shared folders: * ? " < > | ;

When you enter a path that includes an unsupported character, an error message appears. Folder names that include unsupported characters cannot be selected in browse results. You can rename the shared folder to exclude the unsupported character or choose another folder that does not include one of these characters in its name.

Entering Log-in Credentials

When you export an image to a shared folder on a network drive or a stand-alone computer, you must enter log-in credentials. After you connect to a shared folder, the connection remains active until you log out.

Before you export images to a shared folder, obtain the following information:


- The UNC path to the shared folder on the network drive or stand-alone computer.
- Log-in credentials to the shared folder

If you do not have write permission to the shared folder, see your system administrator for write permission.

To export images to a shared folder

1. Verify that the imager is connected via an Ethernet cable to a network outlet or to a stand-alone computer.
2. Do one of the following:
 - In the Gallery, select one or more images to export.
 - In Image View, display an image to export.

3. Tap Send/Save in the bottom toolbar.
4. Select export options.
5. Tap Save to Network and then tap Browse.

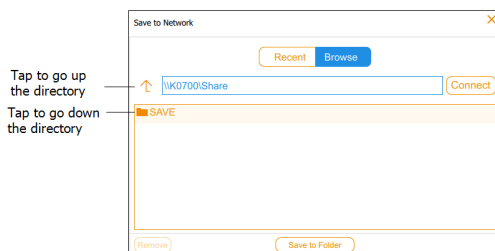

6. Do one of the following:

- Enter the UNC path name to the shared folder on the network drive or stand-alone computer.

Tip: This is the path name you noted when you set up the shared folder. See either [Determining the UNC Path to a Network Folder on a Windows Computer on page 149](#) or [Determining the UNC Path to a Network Folder on a Mac Computer on page 150](#).

- Tap the path name you want under Save in the Save to Network dialog box.

7. Tap Connect.

If you are not connected to the network drive or computer, the system prompts you for log-in credentials.

8. Enter the log-in credentials to the shared folder and then tap OK.

For a network drive, type the network domain name, your user name on the domain, and your domain password in this form:

domain_name\username

for example, **Global.xyz.com\jsmith**

9. Tap Save to Folder. A progress box monitors the image export.
10. When the export finishes, tap OK.


Exporting to a Recent Location

When you export images to a shared folder on a network drive or stand-alone computer, the software saves the location you enter. The next time you export an image, the Save to Network dialog box displays the ten most recent locations you entered, most recent location first.

To export images to a recent location

1. Verify that the imager is connected via an Ethernet cable to a network outlet or to a stand-alone computer.
2. In the Gallery or in Image View, select the images you want to export.
3. Tap Send/Save in the bottom toolbar.
4. Select export options.
5. Tap Save to Network.

The Save to Network dialog box appears.

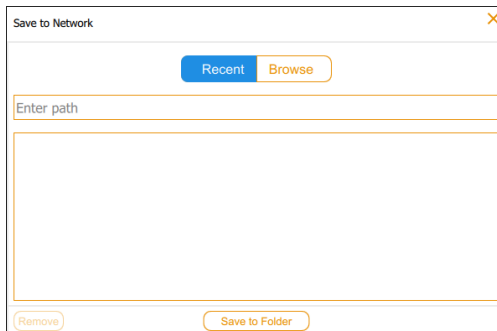The image shows a 'Save to Network' dialog box. At the top left is the title 'Save to Network' and a close button (X) at the top right. Below the title are two buttons: 'Recent' (highlighted in blue) and 'Browse' (highlighted in orange). Under these buttons is a text input field with the placeholder text 'Enter path'. Below the input field is a large, empty rectangular area, likely for a list of recent locations. At the bottom of the dialog box are two buttons: 'Remove' and 'Save to Folder'.

6. Tap a shared folder in the recent locations list and then tap Save to Folder.

A log-in dialog box appears if the connection to the network drive or stand-alone computer is not active.

7. Enter the log-in credentials to the shared folder.

To enter the log-in credentials for a network drive, type the network domain name, your user name on the domain, and your domain password. Type the domain and user name in this form:

domain_name\username

for example; **Global.xyz.com\jsmith**

8. Tap OK.

A progress box monitors the image export.

9. When the export finishes, tap OK.

Removing Network Locations

To remove network locations from the list

- ▶ Tap a location and then tap Remove.

Disconnecting from a Shared Folder

The imaging system remains connected to a shared network folder until you log out or turn off the imaging system or until the network disconnects you.

Chapter 11 System Settings


Setting the System Date and Time

Before you begin to work with the imaging system, check whether the date and time settings are correct for your locale. If your system is connected to a network, date and time are synchronized to the network setting.

The date and time appear in the default names assigned to acquired images. The system date and time are determined by the time zone and the current time settings.

To set the system time zone

1. Tap Settings in the main toolbar.
2. Tap Set Time Zone in the list that appears.

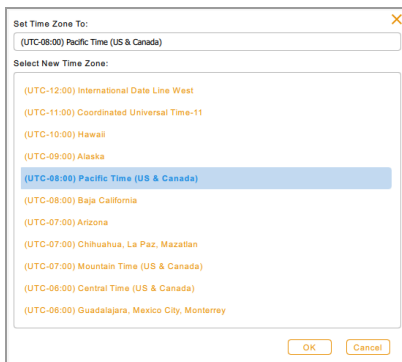


3. In the Select New Time Zone list, swipe to find the time zone for your locale.

4. Tap the time zone to select it and then tap OK.

To set the current Date and Time

1. Tap Settings in the main toolbar.
2. Tap Set Date and Time in the list that appears.
3. In the Calendar box, tap the angle brackets to display the current month.

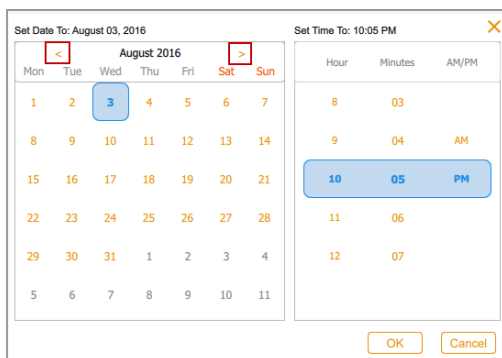

4. Tap the day of the month.
5. Scroll through the Hour and the Minutes lists until the current time appears in the blue band.
6. Scroll to position AM or PM in the blue band under AM/PM.
7. Tap OK.

The selected date and time appear at the top of the all Image Lab Touch screens.

Setting the Sound Volume

You can change the system sound volume or turn off the sound.

To set the sound volume

1. Tap Settings in the main toolbar.
2. Tap Set Sound Volume in the list that appears.
3. Tap a volume level.
4. Tap X to close the list.

Appendix A Maintaining the Imaging System

Reinstalling Image Lab Touch Software

You can reinstall the version of Image Lab Touch currently in use or a previously installed version.

To reinstall Image Lab Touch

1. Obtain the USB flash drive on which the installation software is stored.
2. Verify that the installation software is in the root directory of the flash drive.
3. Ensure that no image acquisitions or exports are in progress.
4. Insert the USB flash drive into the USB port on the imager front panel.
5. Tap Settings on the main toolbar.

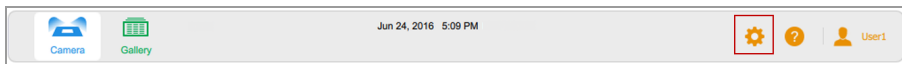


Image Lab Touch detects the installation software in the root directory of the USB flash drive. Update Installation appears in the Settings list.

6. Tap Update Installation and tap OK.


The installation process begins.

Important: Do not interrupt the installation process. When it is complete, the imager restarts automatically.

Maintaining Transilluminator Bulbs

Use 302 nm standard transilluminator bulbs with the imager. (See [Appendix C, Ordering Information](#) for the catalog number.) Depending on the amount of use, the UV bulbs can last for many years.

The six bulbs are organized into three groups of two bulbs each, identified as Bank 1, Bank 2, and Bank 3. When a bulb fails, an error message specifies the bank of bulbs in which a bulb has failed.

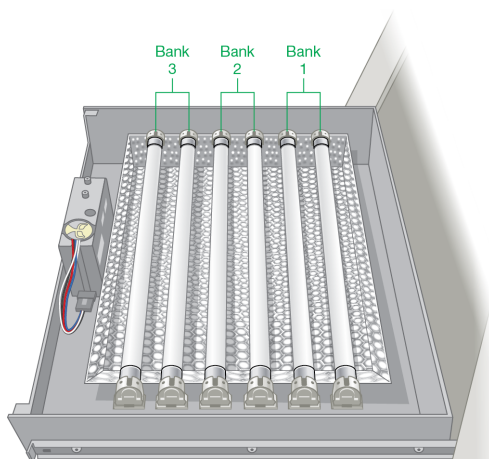

Replacing UV Transilluminator Bulbs

One or more transilluminator bulbs can fail at any time. Bio-Rad recommends that you replace all six bulbs, including those that still work. Replacing all bulbs at once ensures consistent light over the imaging stage.

To replace the transilluminator bulbs, use a 2.5 mm hex wrench.

Important: Transilluminator bulbs contain heavy metals, including mercury. Do not throw used bulbs in the trash. Dispose of them in accordance with local recycling and disposal guidelines.

To prepare to replace the bulbs

1. Turn off the power to the imager.
2. Disconnect the AC power cord from the imager.

To pull out the transilluminator drawer

1. Open the imager door.
2. Use the drawer slide release levers to extend the drawer further and clear the opening. This makes it easier to remove the transilluminator cover in the next procedure.

The levers are located on both sides of the drawer in the drawer slide release mechanism that slides the drawer in and out.

3. Push the lever on the left side down, push the lever on the right side up, and pull.

The drawer extends another centimeter.

To remove the transilluminator cover

1. Loosen the four screws — two on each side of the transilluminator cover.

Tip: Leave the loosened screws in the cover to avoid misplacing them.

2. Remove the transilluminator cover by sliding it forward and then lifting it.


Important: Do not place the transilluminator cover directly on the bench.

3. Place the cover on a nonabrasive surface to avoid scratching or damaging the UV filter glass.

To replace the transilluminator lamps

Important: Use caution when touching the lamps. They can be hot.

1. Rotate a lamp until it loosens and the pins are vertical and aligned with the socket.
2. Remove the lamp.
3. Set the new lamp in place and rotate it until the pins are seated and horizontal.

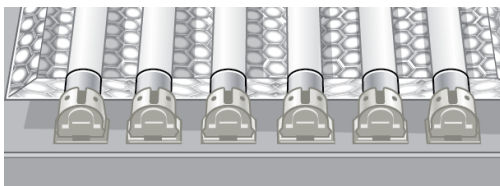

4. Repeat steps 1–3 for each lamp you want to replace.

To replace the transilluminator cover

5. Set the transilluminator cover on the drawer.
6. Secure the cover to the drawer with the screws.

Replacing the Fuses


Replace the fuses with one of the following acceptable replacement fuses:

Name	Number	Voltage	Description
Littelfuse	0217008	250 V, 5 x 20 mm, 8 A	Fast acting
	0217008.HXP		
	0217008.MXP		
Schurter	0034.1425	250 V, 5 x 20 mm, 8 A	Fast acting


To replace the fuses

1. Unplug the power cord from the back of the instrument.


The fuse drawer sits above the power plug.

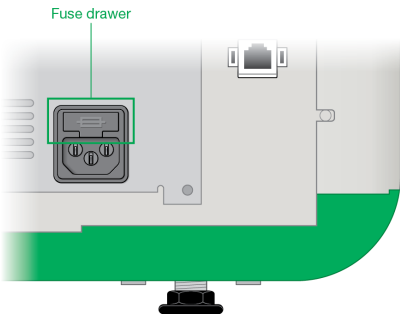


2. Insert the tip of a screwdriver as far as it will go under the center of the fuse drawer.

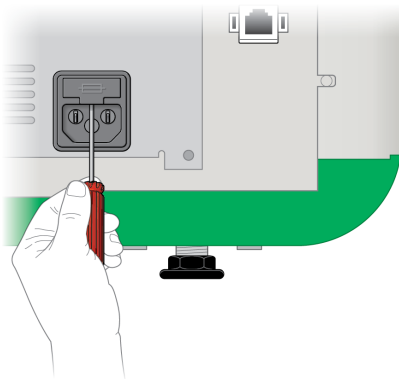

3. Tug hard on the screwdriver to open the fuse drawer.
4. Remove the drawer from the power module.

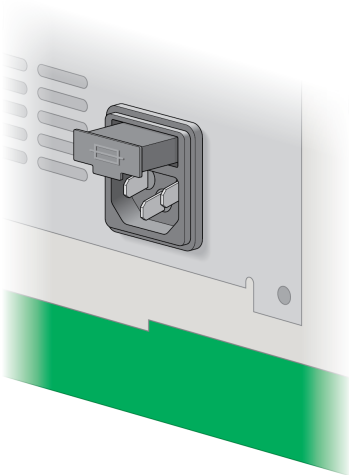

5. Remove the fuses and inspect them to determine whether to replace them.

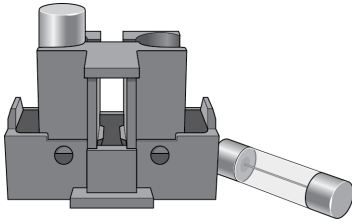

6. Snap new fuses into place.
7. Push in the fuse drawer until it snaps into place.

Appendix B Important Imaging Concepts

This appendix provides additional information on important concepts for working with the imaging systems.

Exposure Settings for Chemiluminescent and Low-Light Applications

You can use the Image Resolution/Time scale with adjustable binning settings to optimize the image you want to acquire. Choose a lower binning setting (higher resolution) for higher-quality images. If bands are faint, choose a higher binning setting for a shorter exposure time and the degree of sensitivity you require.

Image Lab Touch automatically determines the optimal setting required to achieve images suitable for quantitation or publication, as indicated by the Image Resolution/Time Scale. Images acquired at the optimal setting have a resolution of 175 microns or more per pixel. Selecting lower bin settings increases the resolution but might also increase the imaging time. Selecting higher bin settings reduces the resolution and exposure time but might result in a lower resolution than required.

Bio-Rad recommends that you start with the optimal binning setting shown in the Image Resolution/Time Scale, examine the result, and adjust the setting for subsequent images.

Adjusting How Images Are Displayed

Image Lab Touch software optimizes images based on the range of intensity levels in the image and the known behavior of the applications. Use this optimized image as a starting point. Use the histogram scale and the gray scale curve settings in the Transform dialog box to adjust the image brightness and contrast as necessary.

Important: The transform settings change only the *appearance* of the image, not the underlying data.

Note: Chemiluminescent images taken in SAM mode are displayed with a fixed transform so you can compare them. The intensity is set to the bottom third of the full range, and the gamma is set to .75. The fixed transform enables you to distinguish accurately the differences between the images in one SAM acquisition so you can select the best one.

Adjusting Image Brightness and Contrast

Transform adjusts image brightness and contrast, optimizing the image display to highlight features of interest. The minimum to maximum range varies depending on the light and dark values present in the image. Adjustments to brightness and contrast do not change the data. They change only the way the data are displayed.

The frequency distribution histogram shows the total data range in the image and the amount of data at each point in the range. The lightest part of the image is set to the minimum intensity, and the darkest is set to the maximum.

Use the Low and High sliders to narrow the displayed gray scale range.

- The High indicator determines the intensity value shown at the maximum value of the gray scale in the gel image.
- The Low indicator determines the intensity value shown at the minimum value of the gray scale in the gel image.


- The Gamma slider changes the gray scale curve. A value of 1 is linear. A value of <1 redistributes more of the gray scale to the first half of the intensity values. A value of >1 redistributes more of the gray scale to the second half of the intensity values.

A logarithmic histogram can reveal the presence of intensity values that are otherwise obscured. In images with a large background area, the intensity value of most pixels is that of the background. Often, there are too few pixels at the intensity values of the data peaks of interest to make these peaks readily visible in a linear histogram.

A linear histogram can be useful when intensity values are more evenly distributed. Used in combination with the logarithmic histogram, the linear histogram shows more clearly the relative quantity of pixels at each intensity value.

To change the histogram scale

1. Open an image in Image View and tap Transform.
2. Tap the Log Histogram box to choose the logarithmic scale.

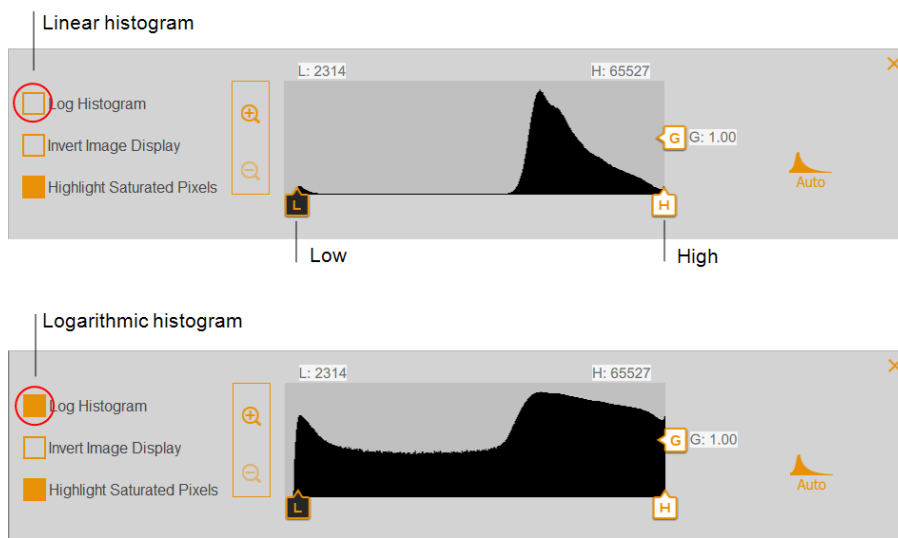


To change the light intensity range displayed in the image

- ▶ Tap and drag the Low or High slider.
 - Drag the Low slider to the right to obscure the background and low-intensity bands and to make high-intensity bands more visible.
 - Drag the High slider to the left to make the background, nonspecific binding, and low-intensity bands more visible. Intense bands might look overexposed.

To adjust the gray scale curve

- ▶ Tap and drag the Gamma slider or touch anywhere in the slider bar.

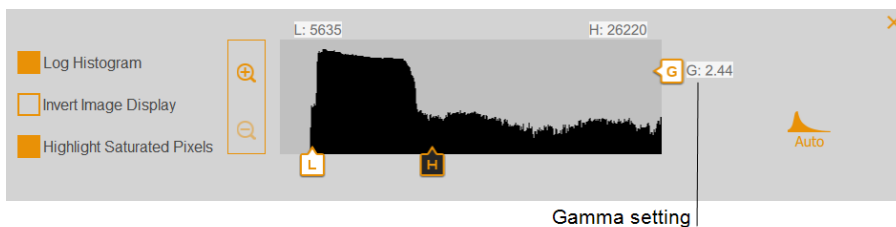

Other Display Options


- **Zoom in and out** — displays greater detail about the intensity range.
- **Highlight saturated pixels** — displays in red areas with saturated signal intensity (higher than a measurable range).
- **Invert image display** — inverts dark and light areas. Clear the box to return to original display.

Restoring Default Settings

After changing the brightness and contrast settings for an image, you can restore the default settings.


To restore default settings

- ▶ Tap Auto.

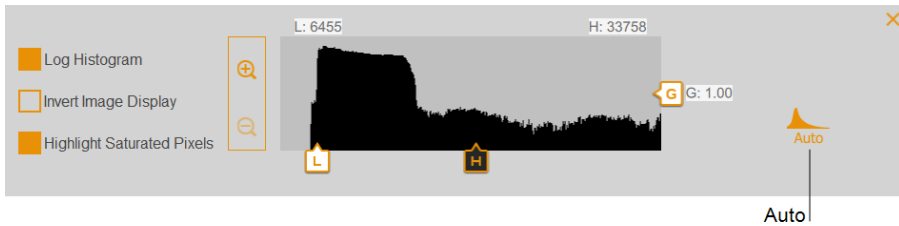




The Image Resolution/Time Scale

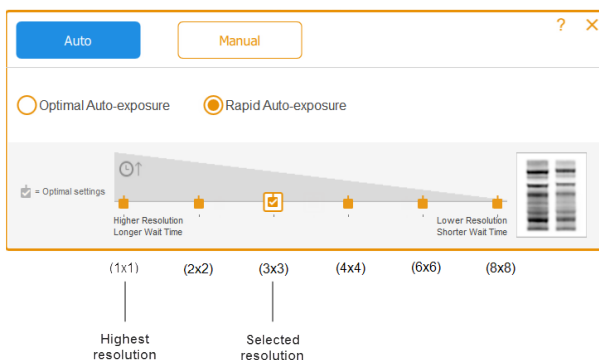The Image Resolution/Time scale includes six pixel resolution settings: 1x1, 2x2, 3x3, 4x4, 6x6, and 8x8. The optimal binning setting required to achieve high-quality images is indicated by a check mark on the Image Resolution/Time scale.




Legend

-  Binning level
-  Selected binning setting
-  Optimal binning setting

Tap a different binning setting on the Image Resolution/Time scale to change the binning level.

When selecting acquisition options, you can view the selected binning setting under Exposure in the imaging settings tile, for example:

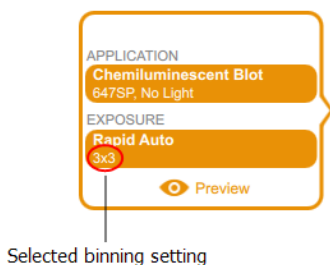

How Binning Settings Affect Image Quality

Pixel binning refers to the process of combining data from adjacent CCD pixels to form a single larger pixel. Doing so results in faster acquisition speeds and, in most images, improved light sensitivity. However, pixel binning is achieved at the cost of reduced resolution. A binning of 2x2 means that 4 adjacent pixels are combined into one larger pixel. A binning of 3x3 means that 9 pixels are combined into one, 4x4 means that 16 are combined into one, and so on.

Pixel binning combines adjacent pixels into a single larger pixel:

Binning option	Combined pixels on the CCD chip
None	
2x2 (4 pixels = 1)	
3x3 (9 pixels = 1)	
4x4 (16 pixels = 1)	
6x6 (36 pixels = 1)	

Using pixel binning, you can balance resolution and exposure time to acquire an image with the sensitivity best suited for your purposes:

- **Sensitivity** — as binning increases, sensitivity to light increases in most images, which improves the ability of the camera to detect faint bands. A binning of 4x4 uses 16 pixels compared to a binning of 2x2, which uses four pixels. As a result, a binning of 4x4 is four times more sensitive than a binning of 2x2.

Note: At higher binning settings, sensitivity might depend on the signal intensity and size of the feature being imaged relative to the background. If the bin contains more pixels than the feature (for example, a band), the intensity is averaged over the binned area. In some images, a higher binning setting might render the feature not visible above the background.

- **Resolution** — as pixel binning increases, fewer pixels appear in the resulting image. As a result, resolution decreases at the higher binning settings.
- **Exposure time** — as pixel binning increases, it takes less time to acquire the image because larger pixels collect more light.


Tip: Zooming in on the area to be imaged increases the resolution and might increase the optimal binning. The loss in resolution with a higher binning setting can be offset by the smaller imaging area and result in a shorter exposure time.

In general, a 2x2 binning setting produces an image that is relatively higher in resolution and requires longer exposure times to achieve equivalent sensitivity to a higher binning setting. An 8x8 binning setting usually produces an image that is relatively lower in resolution and higher in sensitivity (faint bands are optimized), and takes less exposure time to produce.

Tip: For a first acquisition of a sample, you might use a higher binning setting (for example, 8x8) with a shorter exposure time. Evaluate the results and then use a 2x2 binning setting with a longer exposure time on the next acquisition. With some acquisitions the 2x2 binning setting might produce an image with optimal resolution and acceptable sensitivity.

The two gel lanes to the right of the image resolution/time scale show how an image looks given the selected resolution setting. As you tap on the different binning settings, the image displayed changes to show the relationship between the resolution and sensitivity you select.

More intense bands with higher resolution appear at lower binning settings, for example:

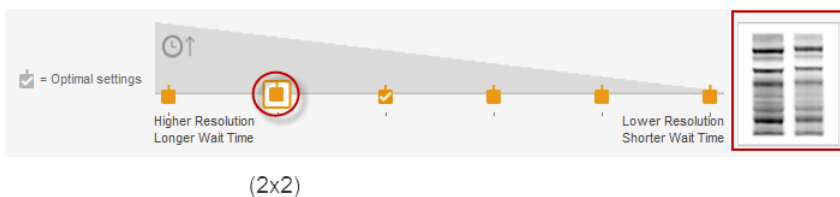

At higher binning settings, faint bands appear more prominent but resolution is reduced, for example:



Appendix C Ordering Information

Catalog #	Description
Imaging Systems	
12003154	ChemiDoc MP Imaging System with Image Lab Touch Software
12003153	ChemiDoc Imaging System with Image Lab Touch Software
Analysis Software	
1709690	Image Lab Software
1709691	Image Lab Software, Security Edition, 1 license
1709692	Image Lab Software, Security Edition, 5 licenses
1709693	Image Lab Software, Security Edition, 10 licenses
ChemiDoc MP/ChemiDoc Imaging System Trays	
12003026	White Tray
12003027	Blue Tray
12003028	Blot/UV/Stain-Free Tray
Optional Accessories	

Catalog #	Description
12003914	UV Safety Shield for ChemiDoc MP/ChemiDoc Imaging Systems
1708377	Holder for Sample Trays and UV Shield
1708185	XcitaBlue Viewing Goggles
12003915	Gel Alignment Template Kit
1708089	Mitsubishi Printer , 100/240 V, USB
1703759	Bio-Rad Fluorescent Ruler
1703760	Gel Cutter Ruler
12005137	ChemiDoc MP/ChemiDoc Imaging System IQ/OQ Kit
Replacement Parts	
1001361	UVB Lamp , 302 nm, 1 each
10026840	Fuse , 8 A, 250 V, 5 x 20 mm, 1 each
1708097	302 nm Lamp Kit , 6 lamps
1707581	Mitsubishi Thermal Printer Paper , 4 rolls
Nucleic Acid Standards	
1708351	EZ Load 20 Base Pair Molecular Ruler
1708352	EZ Load 100 Base Pair Molecular Ruler , 500 µl, 100 applications
1708353	EZ Load 100 Base Pair PCR Molecular Ruler , 500 µl, 100 applications

Catalog #	Description
Protein Standards	
1610373	Precision Plus Protein All Blue Standards , 500 µl, 50 applications
1610363	Precision Plus Protein Unstained Standards , 1 µl, 100 applications
1610376	Precision Plus Protein WesternC Blotting Standards , 250 µl, 50 applications
Buffers	
1610732	10x Tris/Glycine/SDS , 1 L
1610747	4x Laemmli Sample Buffer , 10 ml
Electrophoresis Cells	
1656001	Criterion Cell , includes electrophoresis buffer tank, lid with power cables, 3 sample loading guides
1658004	Mini-PROTEAN Tetra Cell for Mini Precast Gels , 4-gel vertical electrophoresis system, includes electrode assembly, companion running module, tank, lid with power cables, mini cell buffer dam
Blotting System	
1704155	Trans-Blot Turbo Starter System , blotting instrument, includes base, 2 cassettes to hold 1–2 midi or up to 4 mini blotting sandwiches, blot roller, and starter consumable kit
1704156	Trans-Blot Turbo Transfer Pack , mini, PVDF, pkg of 10

Catalog #	Description
1704157	Trans-Blot Turbo Transfer Pack , midi, PVDF, pkg of 10
1704158	Trans-Blot Turbo Transfer Pack , mini, nitrocellulose, pkg of 10
1704159	Trans-Blot Turbo Transfer Pack , midi, nitrocellulose, pkg of 10
1704270	Trans-Blot Turbo RTA Transfer Kit , mini, nitrocellulose
1704271	Trans-Blot Turbo RTA Transfer Kit , midi, nitrocellulose
1704272	Trans-Blot Turbo RTA Transfer Kit , mini, PVDF
1704273	Trans-Blot Turbo RTA Transfer Kit , midi, PVDF
1704274	Trans-Blot Turbo RTA Transfer Kit , mini, LF PVDF
1704275	Trans-Blot Turbo RTA Transfer Kit , midi, LF PVDF
Detection Reagents	
1705060	Clarity ECL Substrate , 200 ml, 25 blots
1705062	Clarity Max ECL Substrate , 100 ml, 15 blots



**Bio-Rad
Laboratories, Inc.**

*Life Science
Group*

Web site *bio-rad.com* **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 43 01 877 89019
Belgium 32 03 710 53 00 **Brazil** 55 11 3065 7550 **Canada** 1 905 364 3435 **China** 86 21 6169 8500
Czech Republic 36 01 459 6192 **Denmark** 45 04 452 10 00 **Finland** 35 08 980 422 00
France 33 01 479 593 00 **Germany** 49 089 3188 4393 **Hong Kong** 852 2789 3300 **Hungary** 36 01 459 6190
India 91 124 4029300 **Israel** 972 03 963 6050 **Italy** 39 02 49486600 **Japan** 81 3 6361 7000
Korea 82 2 3473 4460 **Mexico** 52 555 488 7670 **The Netherlands** 31 0 318 540 666 **New Zealand** 64 9 415 2280
Norway 47 0 233 841 30 **Poland** 36 01 459 6191 **Portugal** 351 21 4727717 **Russia** 7 495 721 14 04
Singapore 65 6415 3188 **South Africa** 36 01 459 6193 **Spain** 34 091 49 06 580 **Sweden** 46 08 555 127 00
Switzerland 41 0617 17 9555 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311
United Arab Emirates 971 4 8187300 **United Kingdom** 44 01923 47 1301

